Service Information

Inhalt

Contents
Contenu
Contenido

Einführung

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Without prior permission in writing by Leica Mikroskopie und Systeme GmbH, this document shall not be reproduced as a whole or in part, by mechanical, photographic, electronic or other means (including conversion into or transmission in machine-readable form); stored in any retrieval system; used for any purpose other than that/those for which it is intended; nor made accessible or communicated in any form to any third party not expressly authorized by Leica Mikroskopie und Systeme GmbH.

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The information in this manual is subject to change without notice.

This manual and its descriptions of functions, data and illustrations are intended to help the user understand the instrument and its use, but imply no claim to accuracy or completeness in detail.

No claim of any kind shall be admissible on the grounds of anything contained in this manual or omitted therefrom.

Exclusion of liability for consequential damage

Leica Mikroskopie und Systeme GmbH hereby expressly excludes the admissibility of any claim for damages or compensation on the grounds of injury, loss, or damage that may arise from the use of the present manual, or from the failure or inability - for whatever reason - to use or understand it properly. This exclusion applies without limitation even if the user or its representative or agent has notified Leica Mikroskopie und Systeme GmbH of the risk or likelihood of such damage, and shall include any claims for damages and/or compensation for loss of profits, operational shutdown, or other financial loss.

Only properly trained, skilled specialist personnel are authorized to serve or repair the open instrument and/or its live components, or perform other work thereon.

The mere possession and/or use of this manual does not convey any qualification upon its user as an authorized, properly trained, skilled specialist. Such qualifications are obtainable only by attendance at and successful completion of service-training courses organized by Leica Mikroskopie und Systeme GmbH. This manual forms part of the basic documentation issued for and used in such courses.

The chapters of the service manual presented in this folder are marked with an **[x]** in the table of contents A 01.

All indications of pages consist of a capital letter stating the chapter according to the table of contents A 01 and a number.

The pages are numbered as follows (e.g. 1.1 or 1.2): the number before the dot is the number of the page. Pages with the number 1 behind the dot deal with the basic state of the instrument, component group or procedure concerned. In case of constructive or production alterations a supplementary page will be prepared which will have the next following number behind the dot.

The alterations have been marked on the illustration pages by providing the part number with one or two little lines:

One line underneath the number:

25 Part has been modified,
can be installed into old instruments.

One line on top of the number:

25 Part has been modified,
cannot be installed into old instruments.
(see chapter "Modification instructions")

One line each underneath and on top of the number:

25 Part has been added.
can be also installed into old instruments.
(see chapter "Modification instructions")

One line each beside the number:

[25] Part has been added,
cannot be installed into old instruments.

An individual parts list is belonging to every illustrated page. On the illustrated page the spare parts are marked with numbers, which correspond to those listed in the 1. column of the parts list. Numbers in brackets mark the component group. Numbers in a circle mark a superior component group (instrument or sales article).

Electrical components are marked with a code according to the circuit diagram in the column "Remarks".

The spare parts list includes the part numbers and part names needed in case of a spare parts order and the required quantity per illustrated page.

If several part numbers are listed underneath an illustration number, pay attention to the differences in the columns "Description" or "Remarks".

A little star * behind a number listed in the 1. column of the parts list means that this part is not presented on the illustrated page or that it is not marked with a number.

In most cases you will find the encoded information on the quality of the standard parts' surface in the 7. to 9. number of the part number. When the colours of the instruments or components are different a corresponding surface quality of the parts to be delivered cannot be guaranteed.

In the column "Remarks" letters and numbers give information on:

Component groups
Membership of component groups
Manufacturing states
Possibilities of replacement
Availability
Orders

The letters have got the following meaning:

- a = from serial number ...
= from state of manufacture ...
= from date/year of manufacture ...
- b = up to serial number ...
= up to state of manufacture ...
= up to date/year of manufacture ...
- E = Replacement for current no. ...
- f = is required for current no. ...
- G = Component group
= this article consists of several parts.
The next following numbers indicate the items (current no. or page no. and current no. of the illustrated page) belonging to this component group.
- i = included in current no. ...
- k = to complete this item current number (s) ...is (are) ...required
- m = to be assembled with current no. ...
- N = no longer available
- NE = no longer available, replacement see current no. ...
- o = but without current no. ...
= this part number does not include the items...
- S = part with the serial number engraved
= part with special engraving
Ordering these parts the serial number or special engraving must be indicated.
The defective engraved part must be enclosed with the order.

- U = the current number(s) is (are) to be replaced with this part
- V = Sales article
The 6-digit part number must be used as long as it is included in the "List of outfits"
- w = by choice - as required
- \$ = electr. component, can be found on the illustrated page underneath the electrical code

The Service Manual describes and illustrates assembly, adjustment, repair and modification procedures. The listed tools can be obtained from the Technical Service.

Those code numbers with an oblique stroke in between are no part numbers. They only serve to help in finding the said part within the spare parts list.

Please note that
before the stroke / behind the stroke
the number of page. the number of illustration.

for example B 1.1/5

is indicated. This example means that the component concerned can be found on page B 1.1, current no. 5.

Individual parts and sub-groups have to be cleaned with unleaded petrol, if possible outside the housing. Lacquered parts and such ones with coloured engraving mustn't be cleaned wet, but only with a dry smooth leather or a brush. Dirty external surfaces with structured enamel finished may be cleaned with an oily rag and rubbed dry.

Parts made of glass and rubber mustn't come into contact with petrol, oil or lubricants.

Such points to be provided with lubricants or adhesive are marked with numbers (311) or letters. These auxiliaries are obtainable from the Technical Service. Any other auxiliaries are not allowed as they do not guarantee a faultless function. The parts to be lubricated must be clean and dry.

Please note that a lot of lubricants do not harmonize with several plastic materials that is to say that they possibly dissolve them or cause their swelling. The lubricants should come into contact with the marked surfaces only. Components to be provided with adhesives have to be free of lubricant. Residual adhesive has to be removed with thinner.

Spare part order

Spare parts have to be ordered by the customer's number and the kind of shipment you prefer. Orders are accepted at department TS-Logistik (Fax no ++49 6441 29 2853). You will find the names and part numbers of the part to be ordered in the spare parts list.

In case of orders for articles that are delivered by the metre, 1 piece means 1 metre.

Let us have a special notice if you do not want a part-delivery. In case of any checkbacks with regard to spare parts orders please refer to the TS-order number and the date of the order concerned.

Sales articles (see "List of outfits") must be ordered only for exchange with the Technical Service Division.

We are not obliged to supply you with all parts and component groups included in the spare parts list. In case of orders for parts that are not available, we reserve the right to deliver such parts that include or replace the one you have ordered. In this case the label on the part and the invoice and/or listing will inform you accordingly.

Safety hints

Before you open the instrument pull off the mains plug and/or disconnect the unit from the mains. If a repair, an adjustment or a maintenance procedure of the open instrument is unavoidable these procedure should only be performed by experts who are familiar with possible dangers.

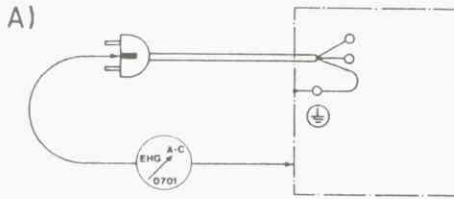
If a harmless operation of the instrument can no longer be ensured, it has to be switched off and secured against unintended starting.

As a matter of principle the national and international safety rules have to be observed.


Remark: Even when the instrument has already been switched off, some internal parts, such as capacitors, might be still current-carrying.

Safety hints


- 1.) After the repair and/or maintenance the instrument must be in accordance with those safety regulations that were in force at the time of the instrument's delivery.
- 2.) In case of maintenance that could be possibly performed by the customer himself, please observe the hints given in the operating instructions.
- 3.) All electronic equipment is subject to safety regulations in accordance with IEC 1010-1 / EN 61010-1 (Safety regulations for electronic surveying and controlling instruments).
Lamps with high-pressure bulbs are subject to IEC 598-1 / EN 60598-1.
- 4.) As a matter of principle all parts, such as transformers, anti-interference filters, which do have direct contact with mains voltage have to be checked after the repair of the instrument and/or exchange of parts.
 - A) Check of the earth conductor connection $\leq 0,3 \Omega$ (with mains cable)
 - B) Measurement of the insulating resistance $R_{isol} \geq 10 M\Omega$
 - C) Measurement of the leakage current:
 $I_{max} = 3,5 mA$ fortified instruments
 $I_{max} = 0,75 mA$ movable instruments
- 5.) For these checks use appropriate testing tools such as EHG-tester VDE 0701.
- 6.) Whenever you carry out leakage current measurement by means of EHG-Tester in accordance with VDE 0701 (B.)
 - a) the current will be higher than in case of measured as test value B, which is caused by the paralleling of the two feeders. Therefore the maximum values for every instrument have to be stipulated.During the measurement the instrument to be tested has to be switched on. Mains connectors have to be bridged.
- 7.) In case of repairs inside circuits (such as the exchange of IC's) carrying signal- or low voltage = 42V, the afore-said tests are unnecessary if all safety regulations have been observed.
- 8.) Anyway, correctly connect the earth conductor connections (green/yellow) and/or put toothed washers underneath the screws when you screw up the parts. If you are not sure, check by means of an appropriate testing instrument!
Moreover do not forget to re-establish all potential compensating lines.
- 9.) Caution !
To prevent malfunctions and/or electrical interference, systems installed or assembled by the user/operator must not exceed the limits specified for class B equipment in the DIN 57871 and VDE 0871 standards.

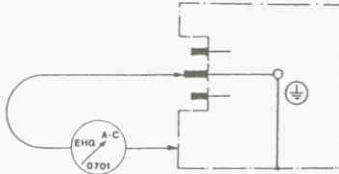
Measurement of the earthed conductor-resistance of instruments with the mains connection cable fixed ($\leq 0.3 \Omega$).

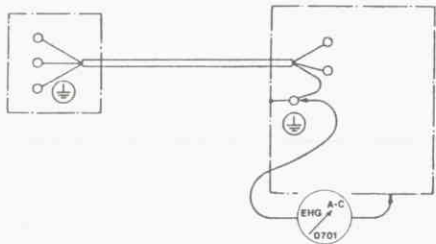Measurement of the earthed conductor-resistance of instruments with detachable mains connection cable ($\leq 0.2 \Omega$).

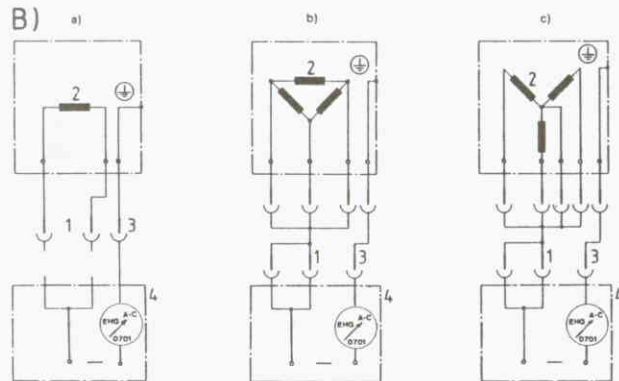
Measurement of the earthed conductor resistance of instruments permanently connected ($\leq 0.2 \Omega$).


- 1 Connection between surveying instrument and the instrument to be checked.
- 2 Internal circuit.
- 3 Connection between surveying instrument and the earthed conductor connection and/or the touchable metal parts
- 4 EHG instrument

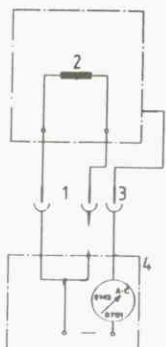Measurement of the insulating resistance; protection category I.

- 1 Connection between surveying instrument and the instrument to be checked.
- 2 Internal circuit.
- 3 Connection between surveying instrument and touchable metal parts.
- 4 EHG-instrument.

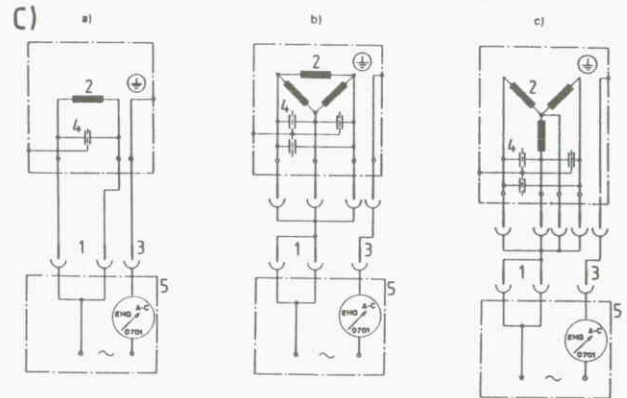Measurement of the insulating resistance; protection category II and III.

- 1 Connection between surveying instrument and the instrument to be checked.
- 2 Internal circuit.
- 3 Connection between surveying instrument and the earthed conductor connection and/or the touchable metal parts.
- 4 Capacitor
- 5 EHG-instrument

Measurement of the leakage-current

How to prevent components from being destroyed by electrostatic charge:

In case of repair please observe the following instructions.


- 1.) Working with components or component groups sensitive to electrostatic charge is only advisable at a workshop place, particularly intended for that purpose. It must be equipped with

- a) a grounded, conductive table coating.
- b) an armlet connected to the table coating.

The pre-resistance of the connection cable between table and ground should be 200kΩ to 1MΩ.

- 2.) Before you start working switch off the instrument and/or pull off the mains plug.
- 3.) Do only insert and/or solder IC's with the armlet put on and the soldering iron grounded.
- 4.) Exclusively use conductive, antistatic special packings with clear markings to pack up spare parts and components after repair and check.

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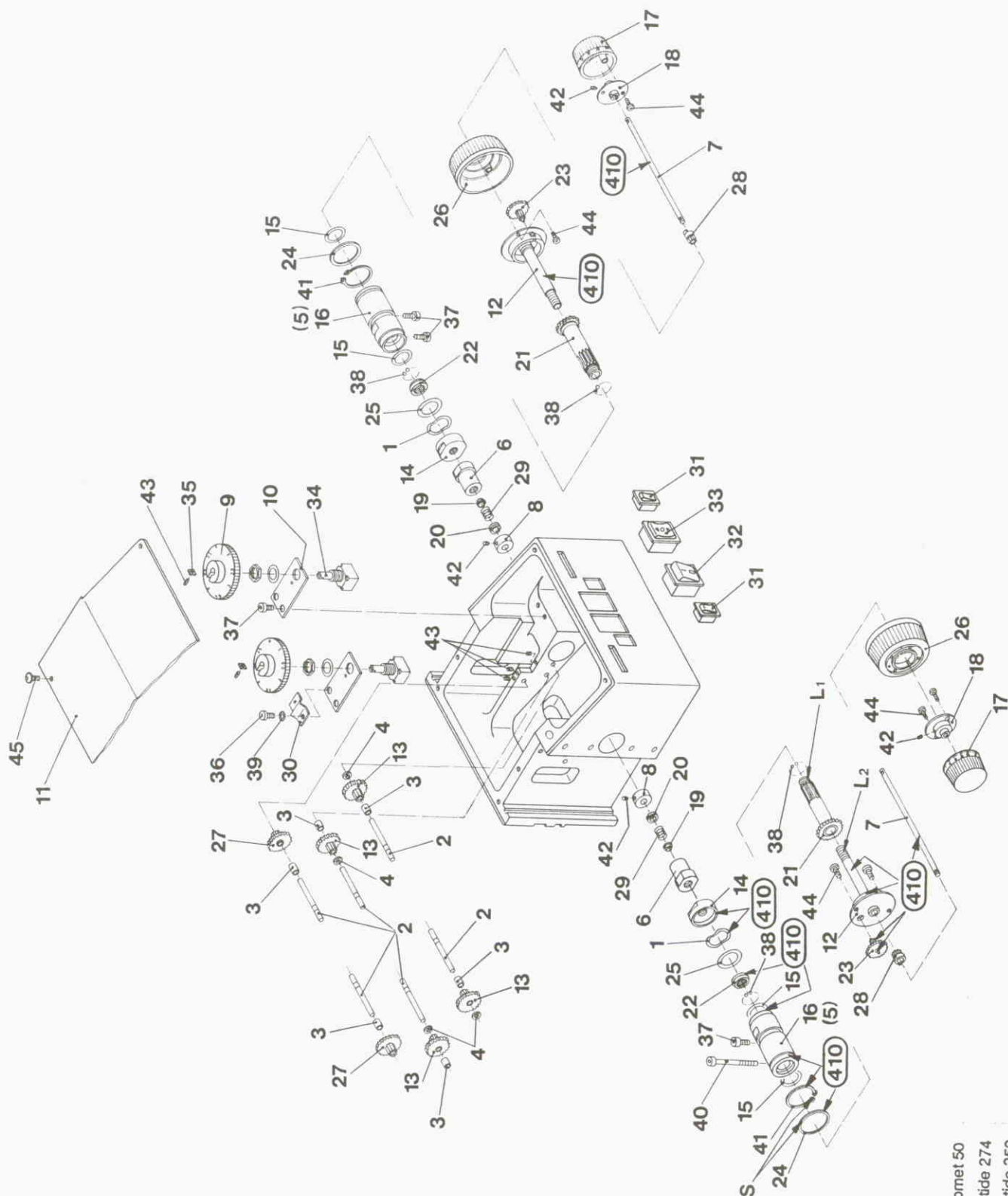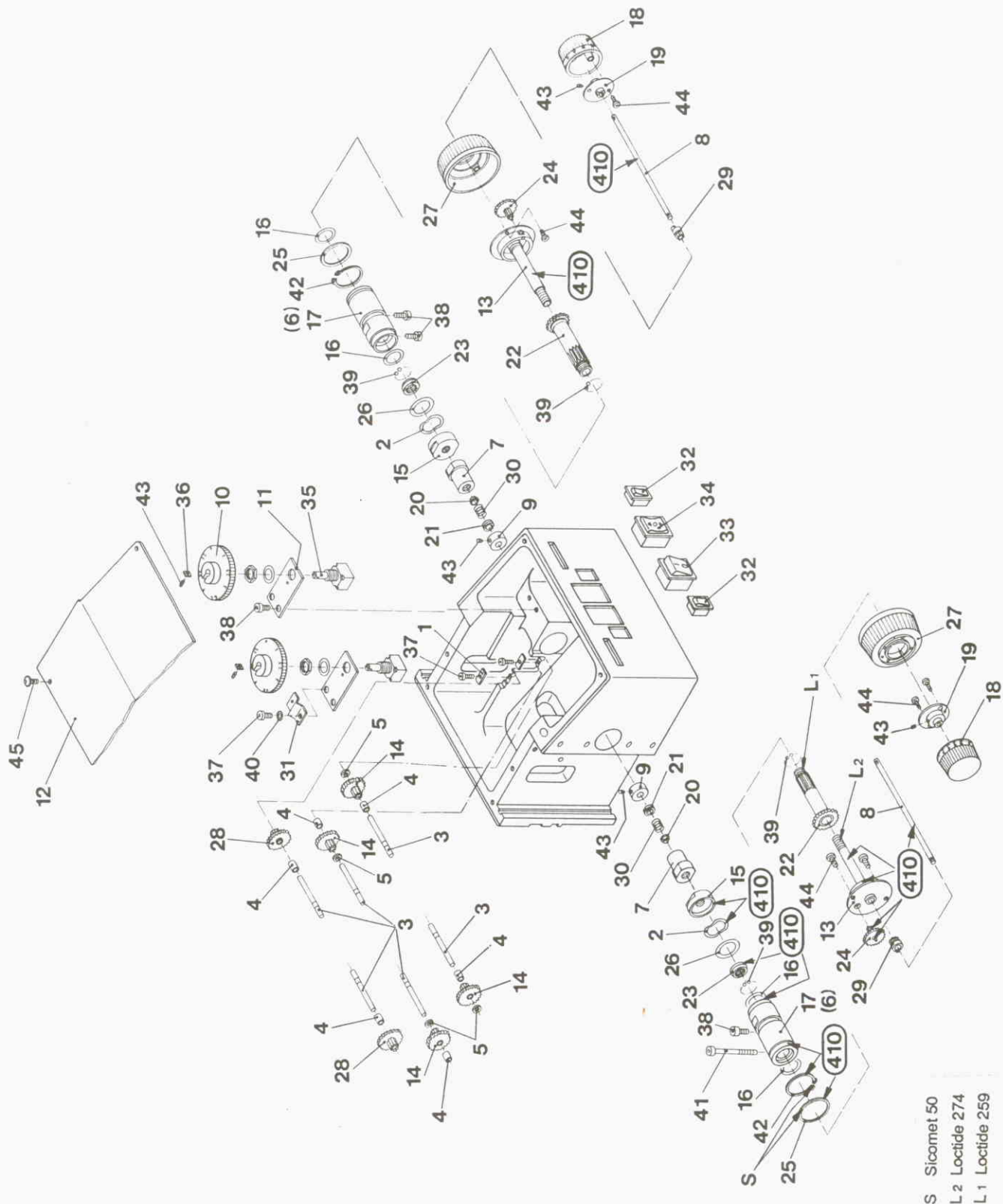
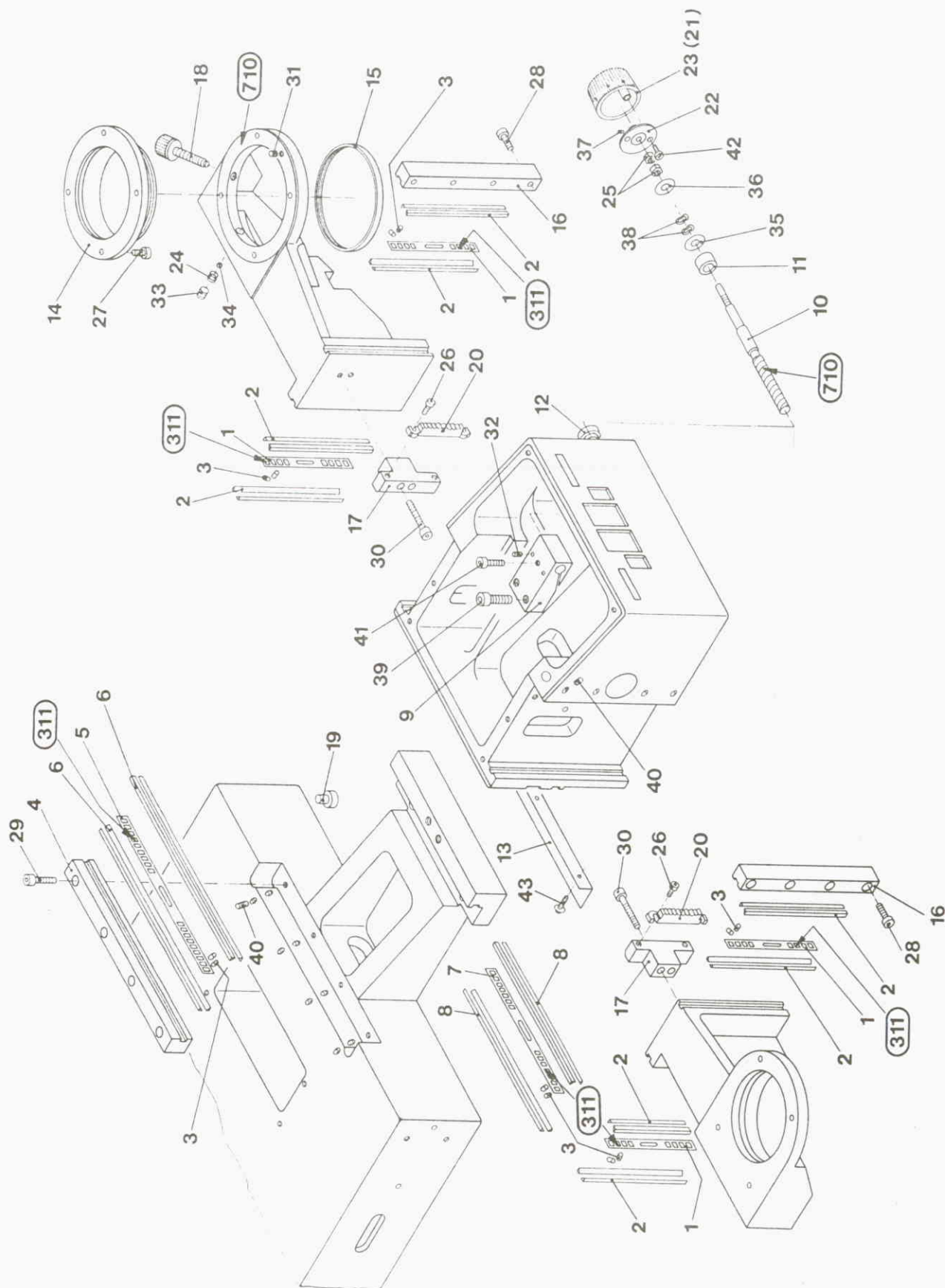
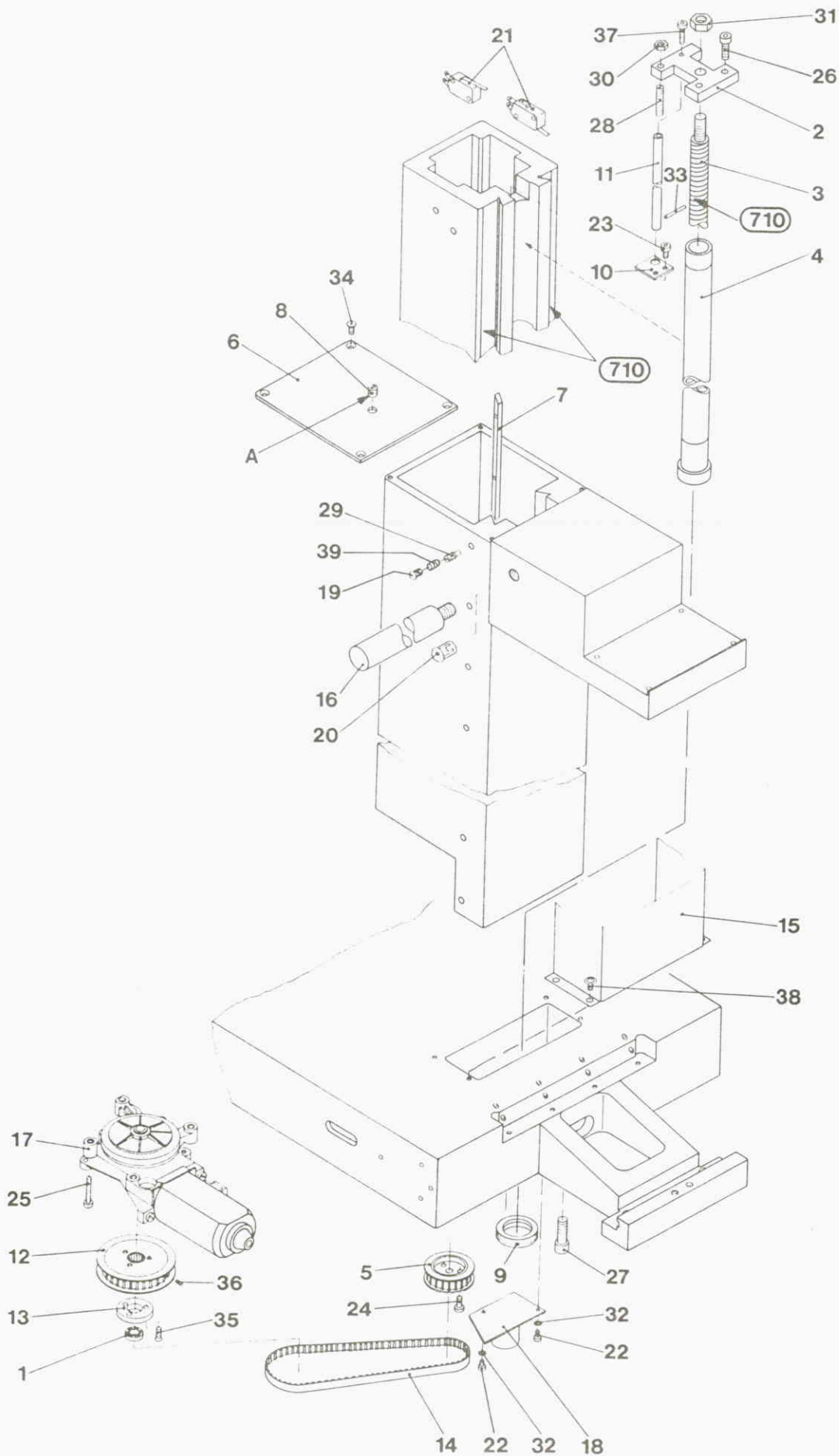
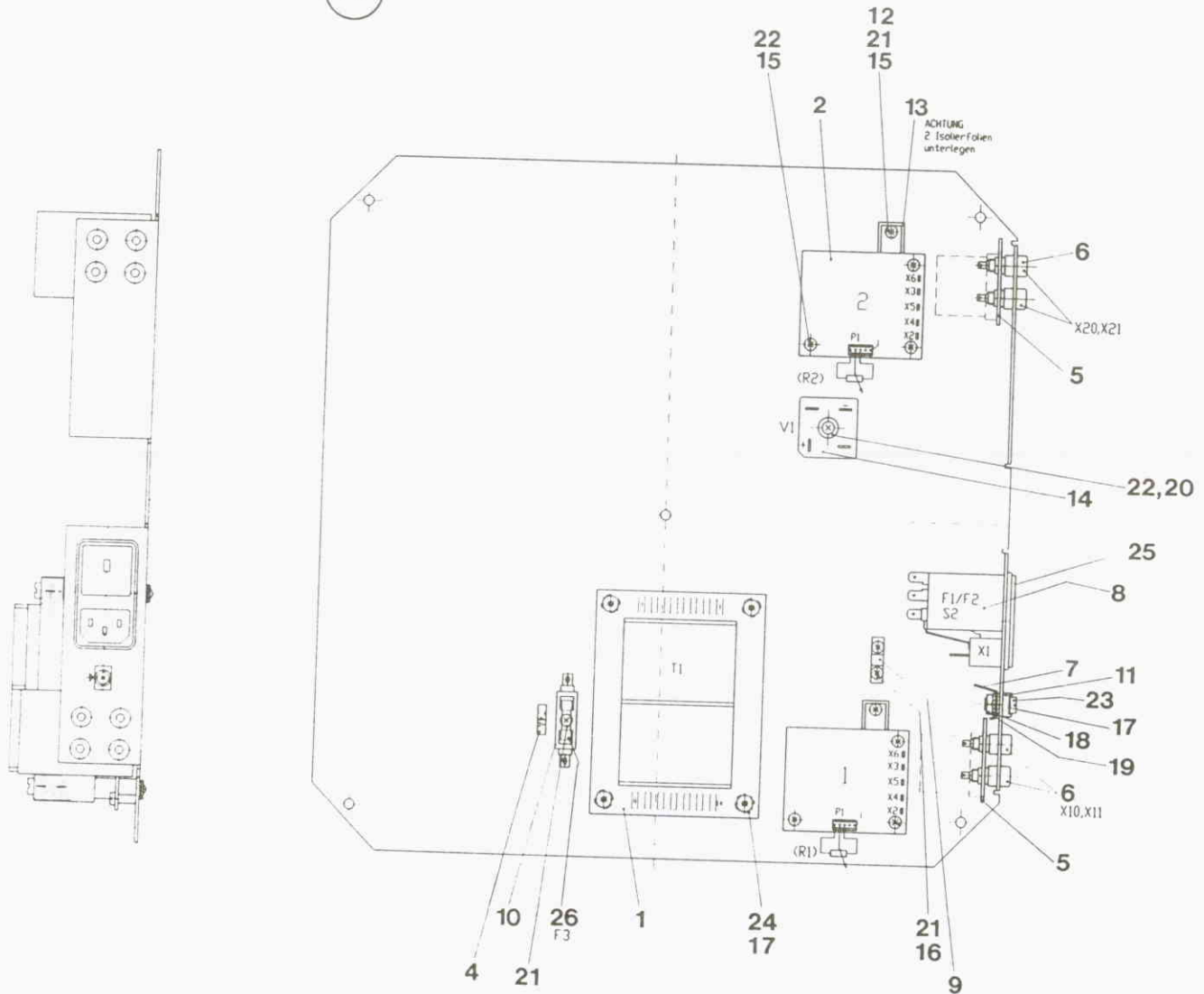
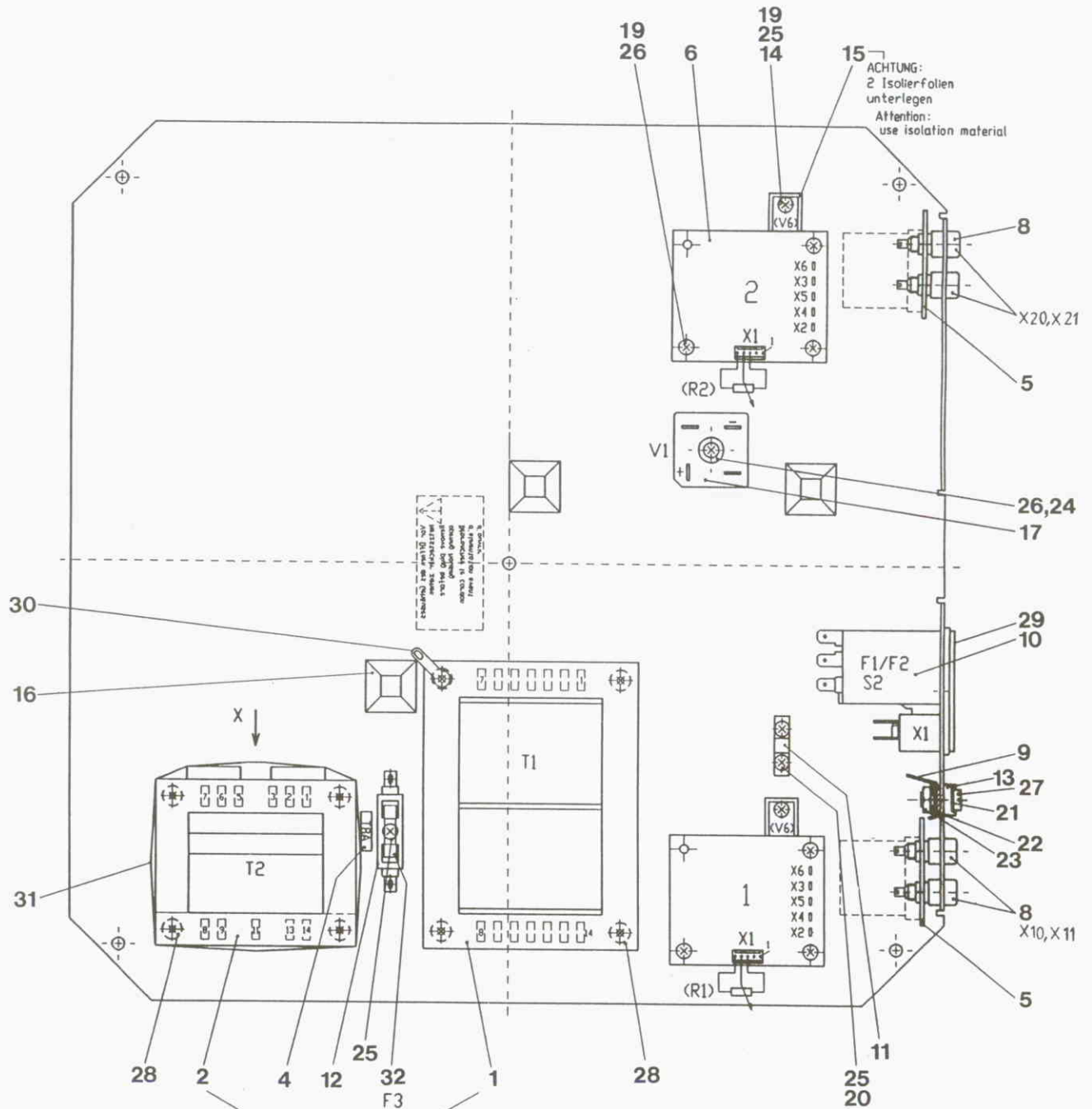
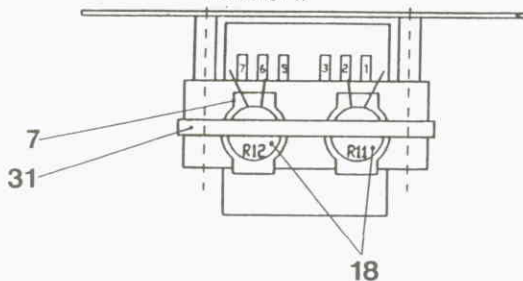
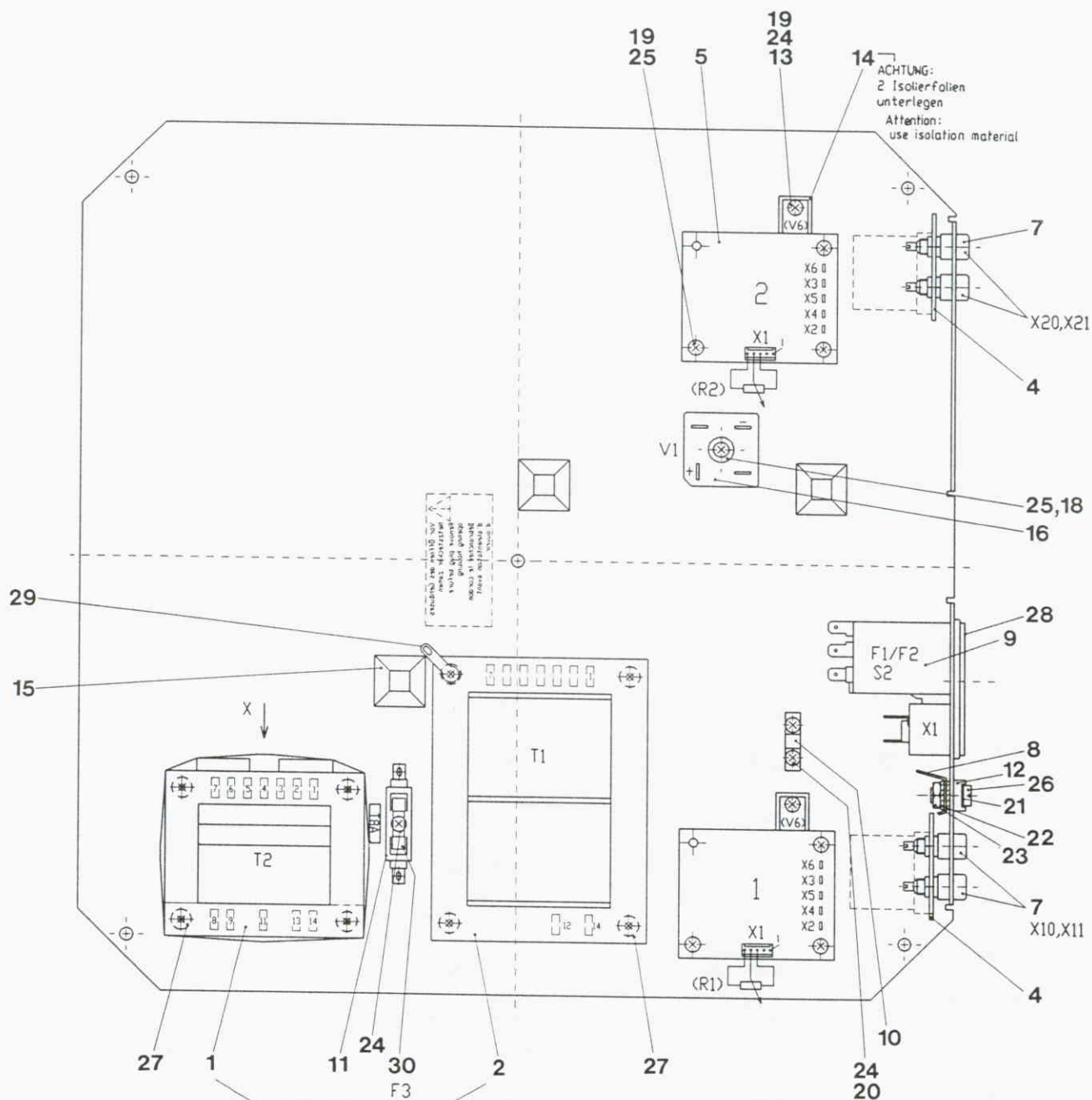
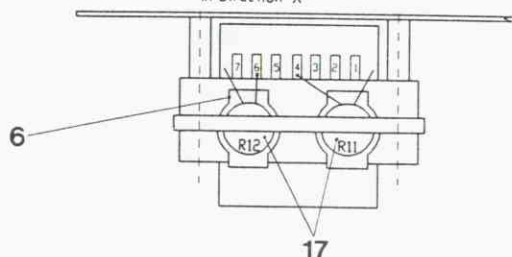
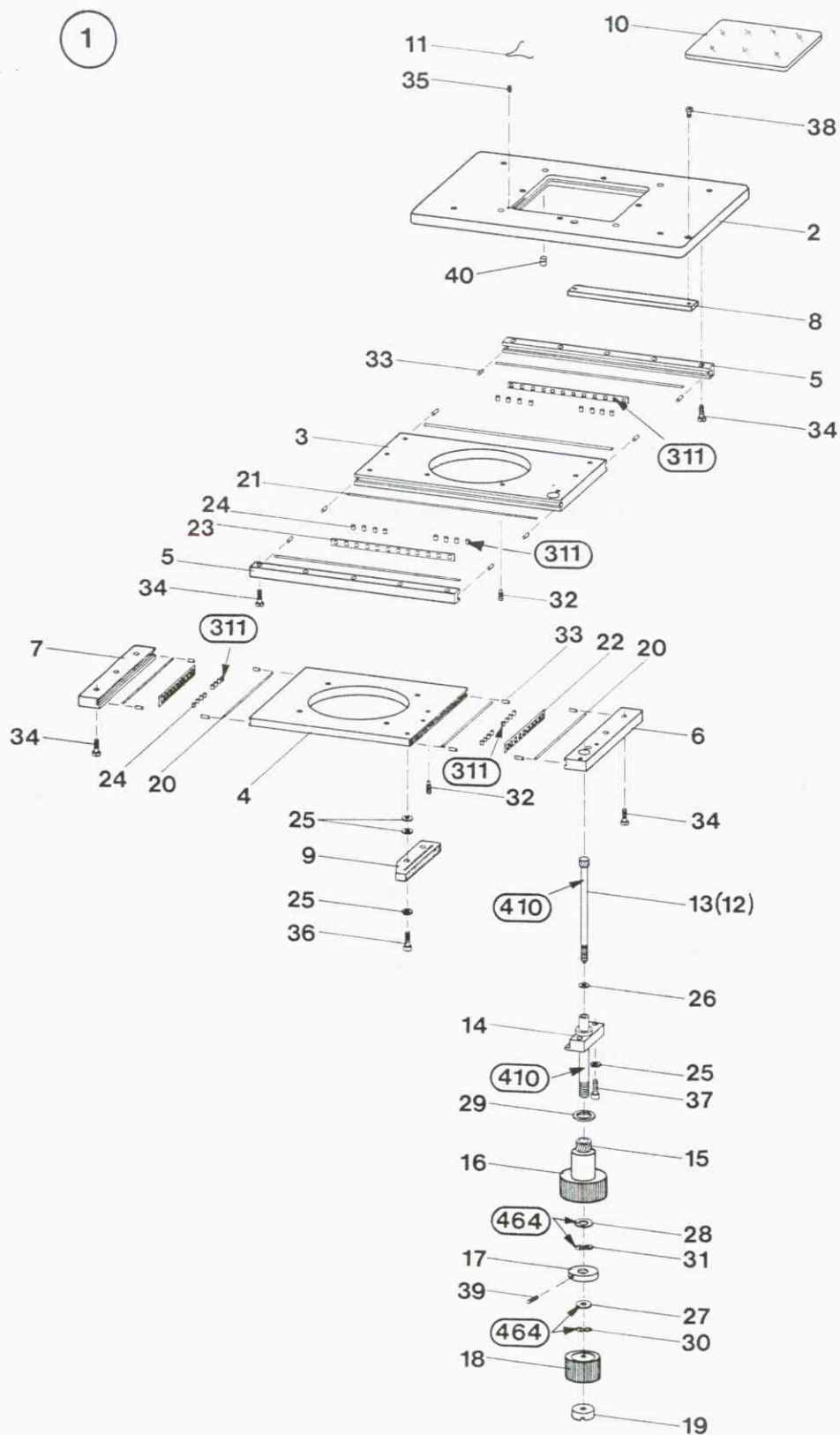
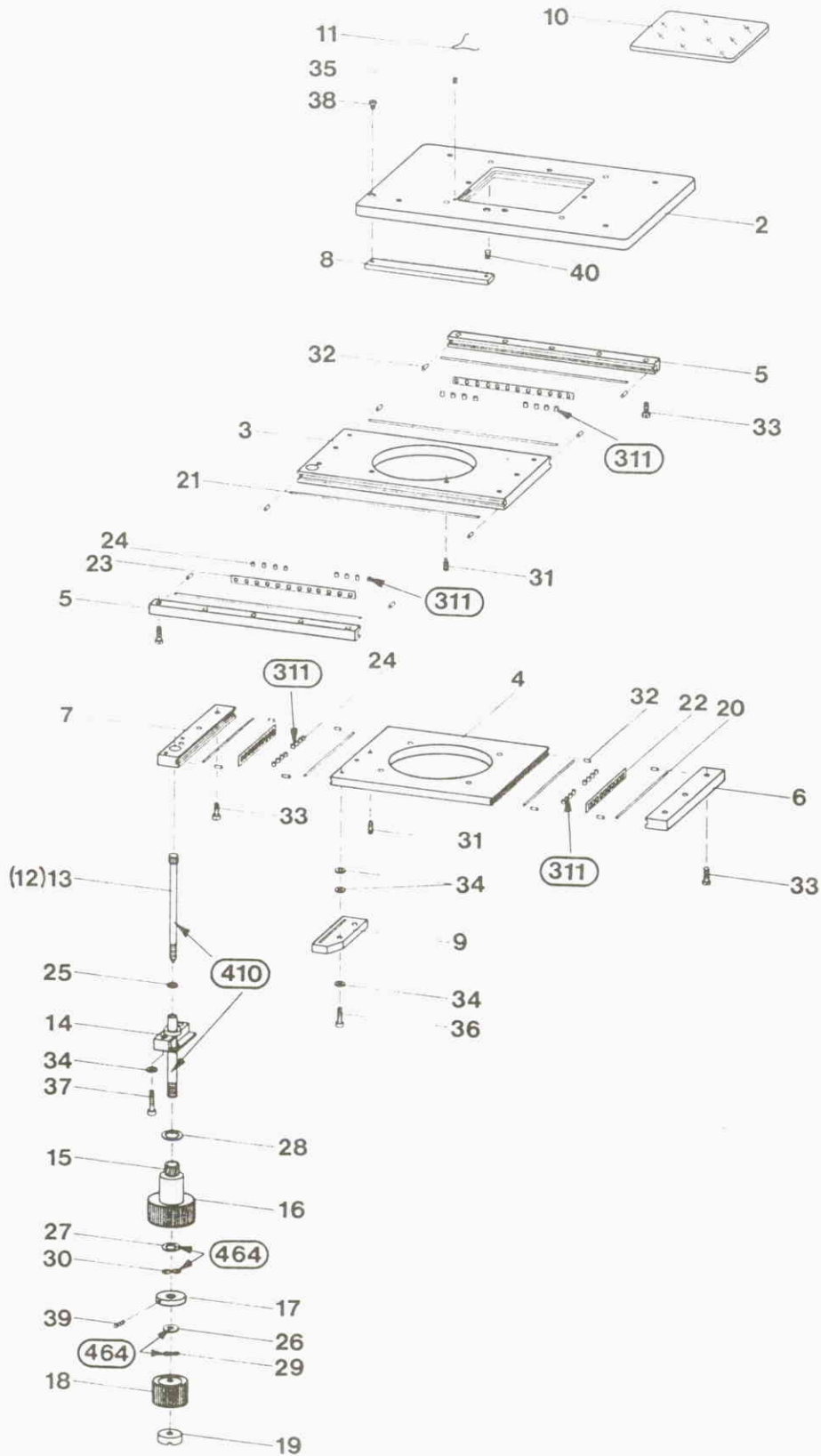
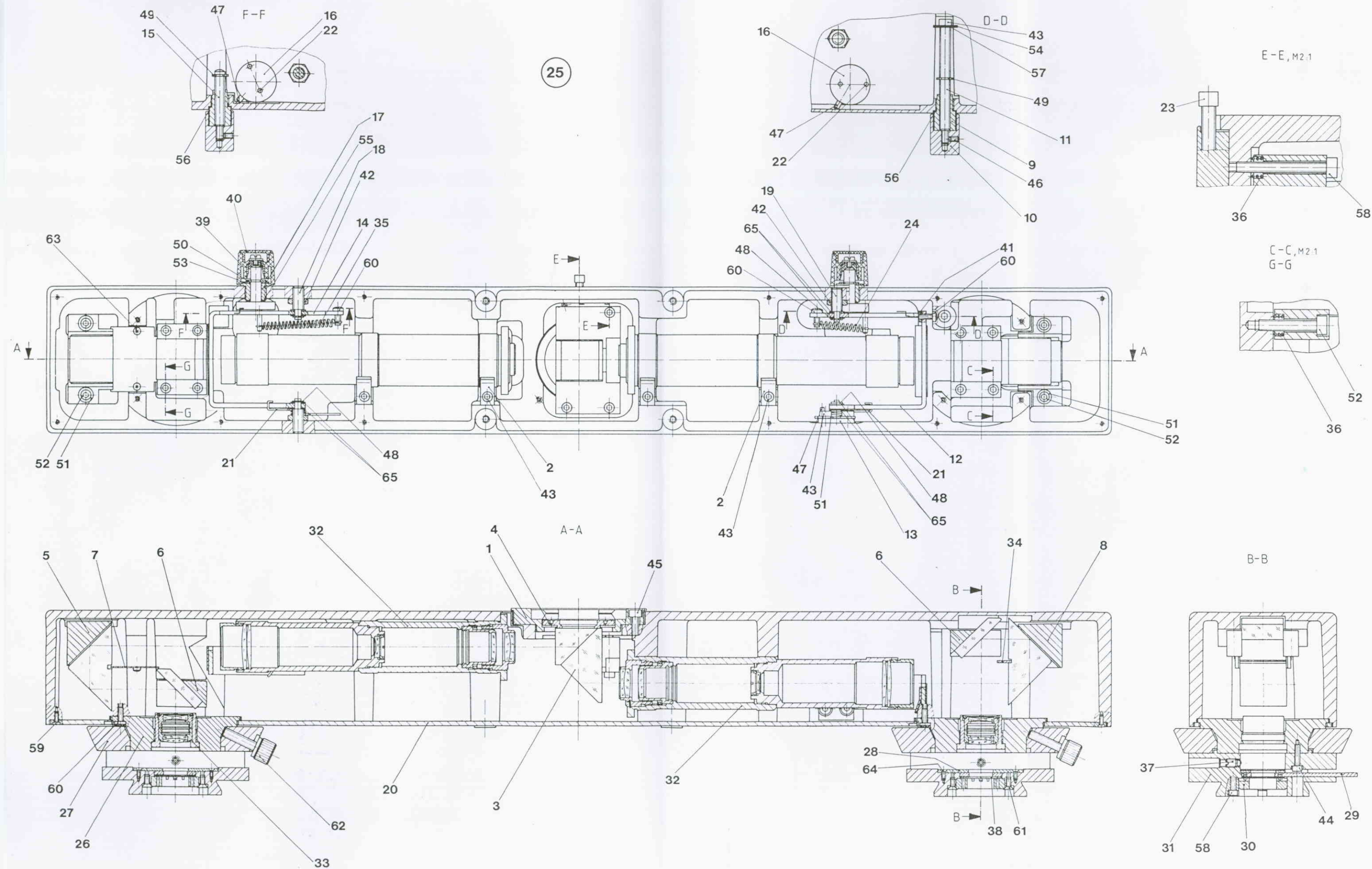
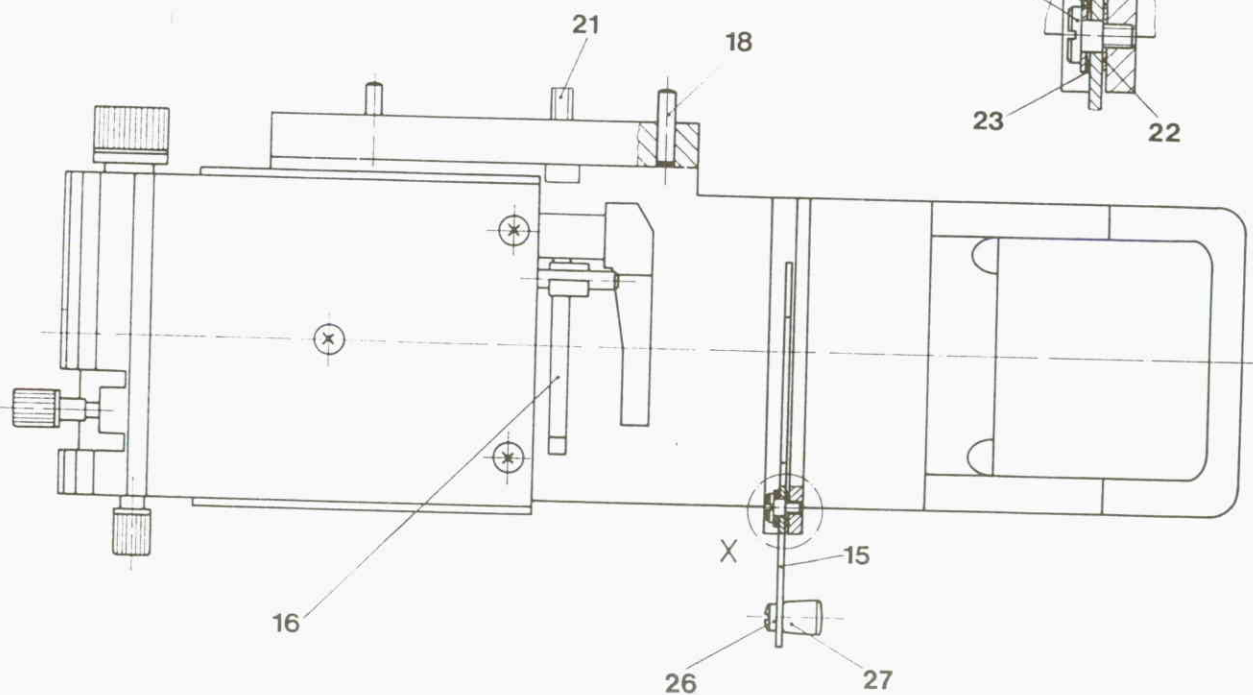
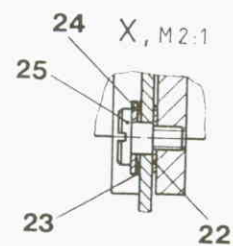
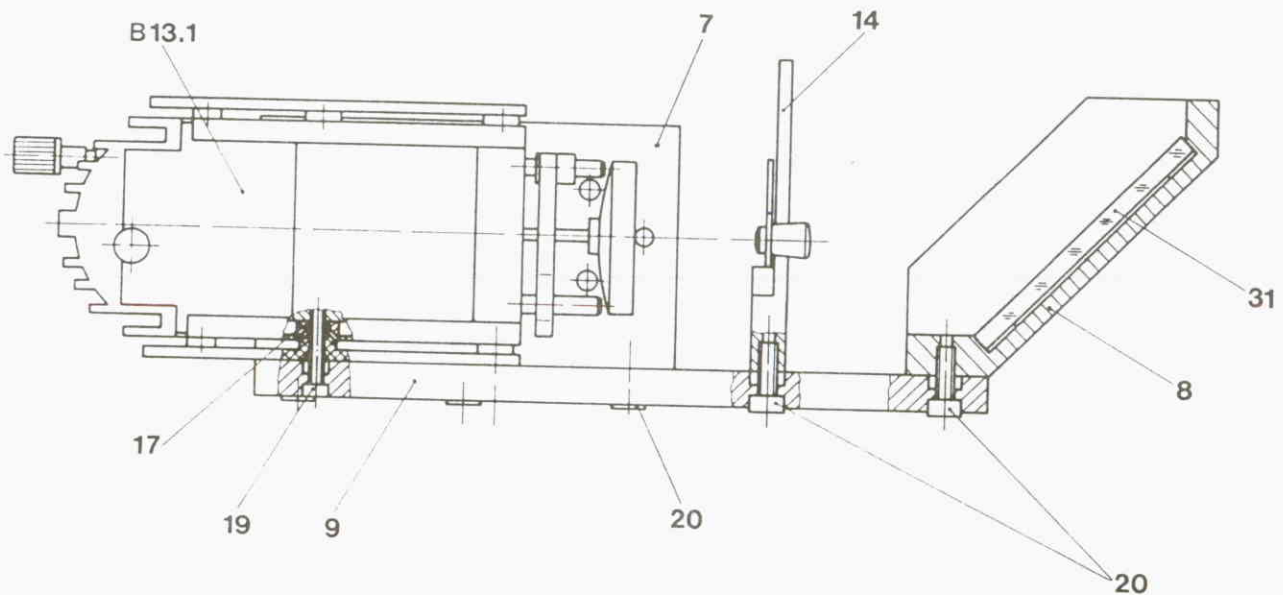
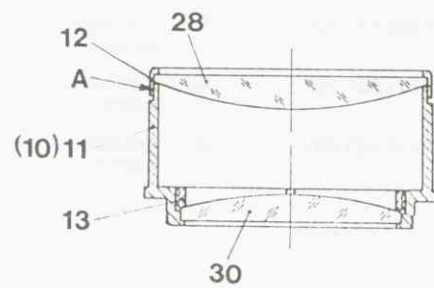
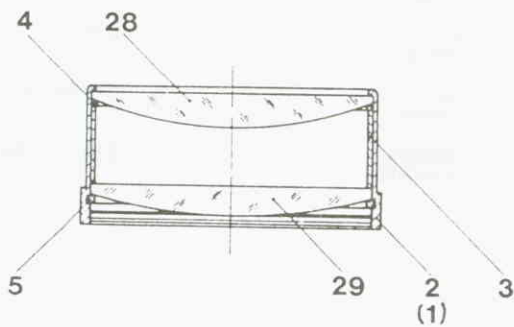
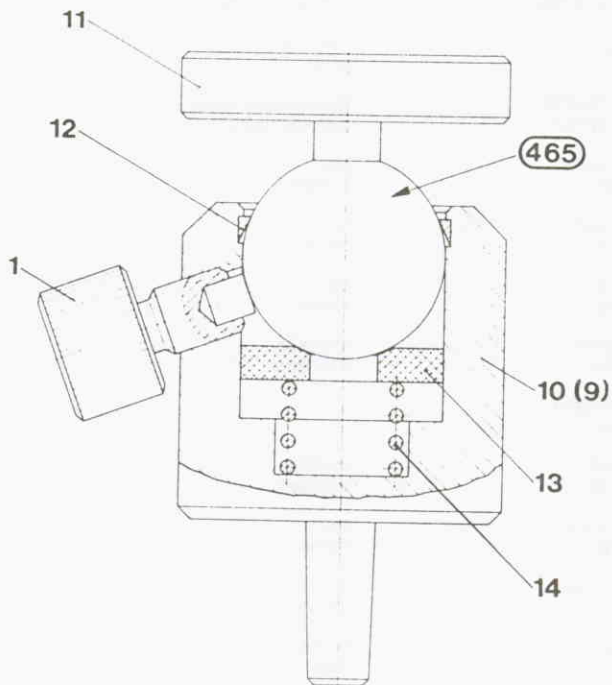
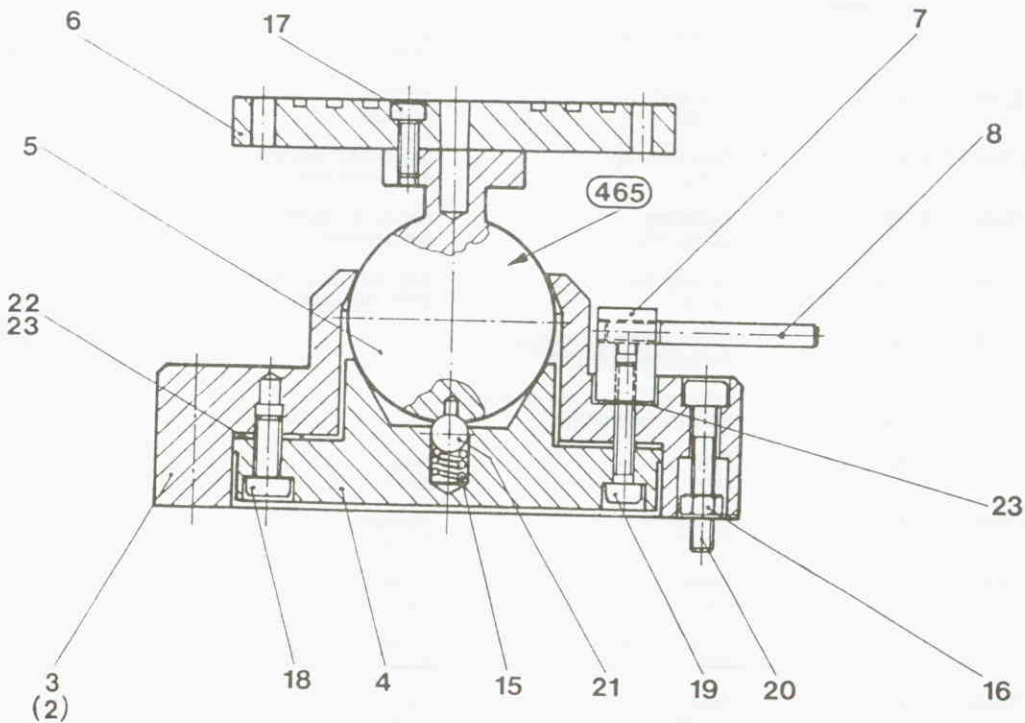
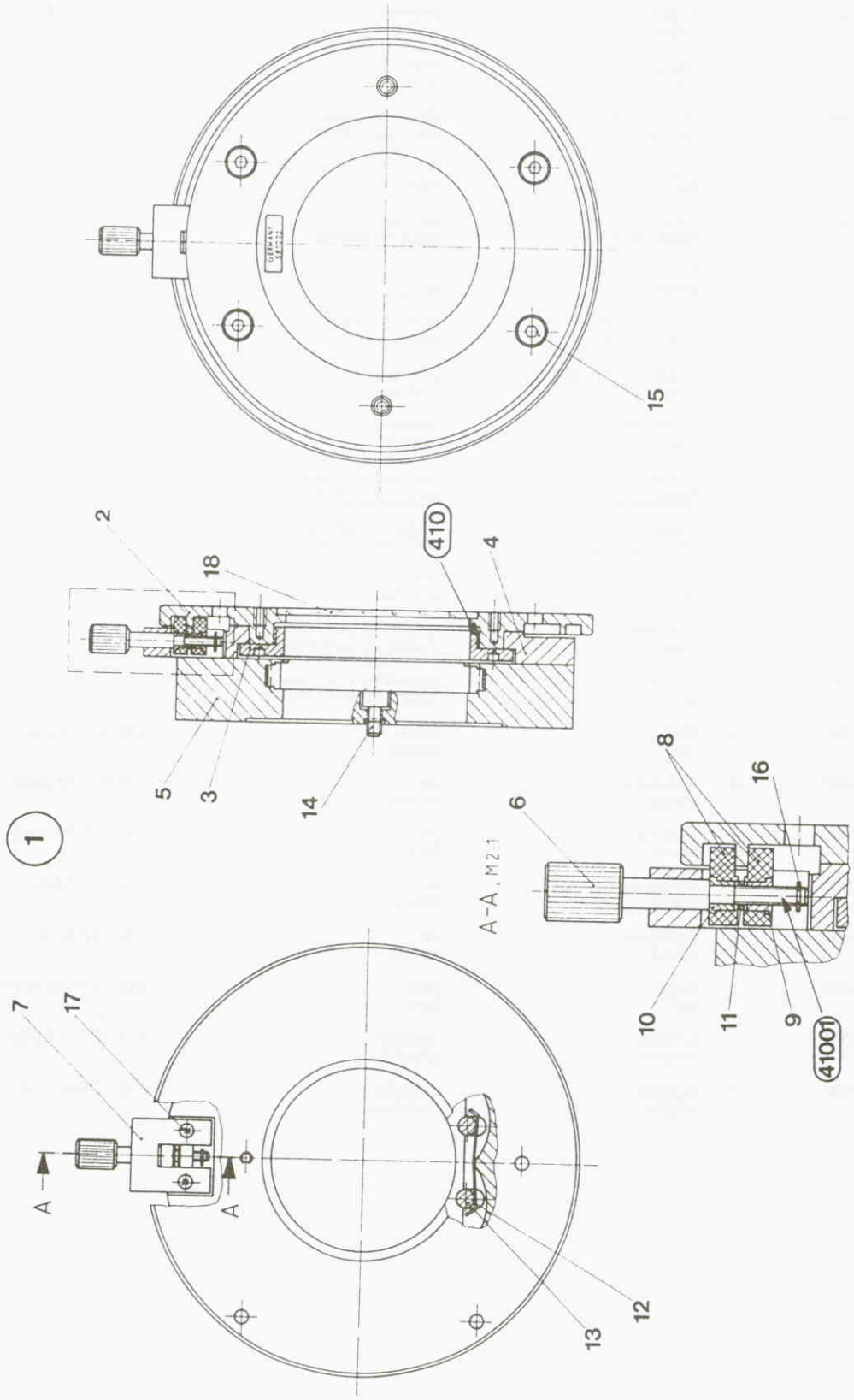
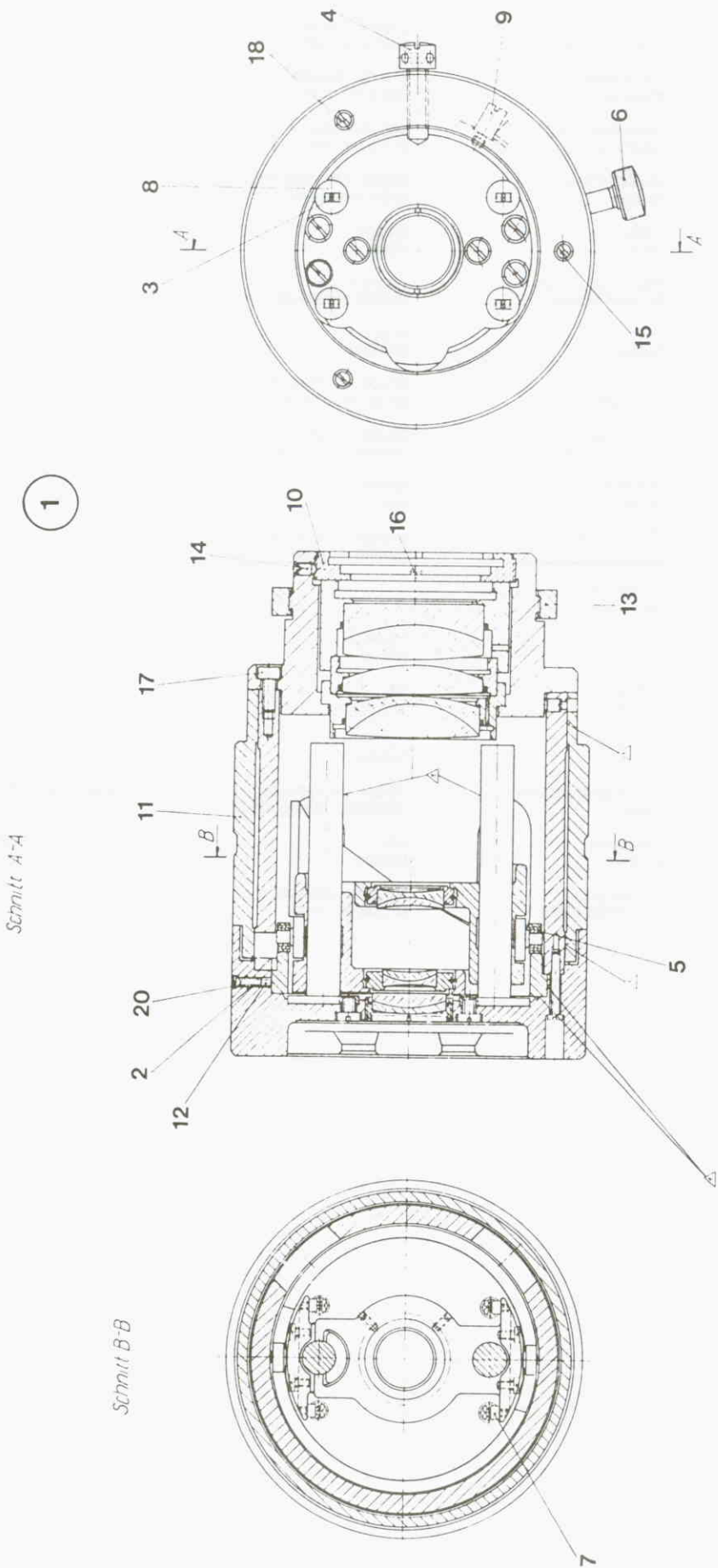
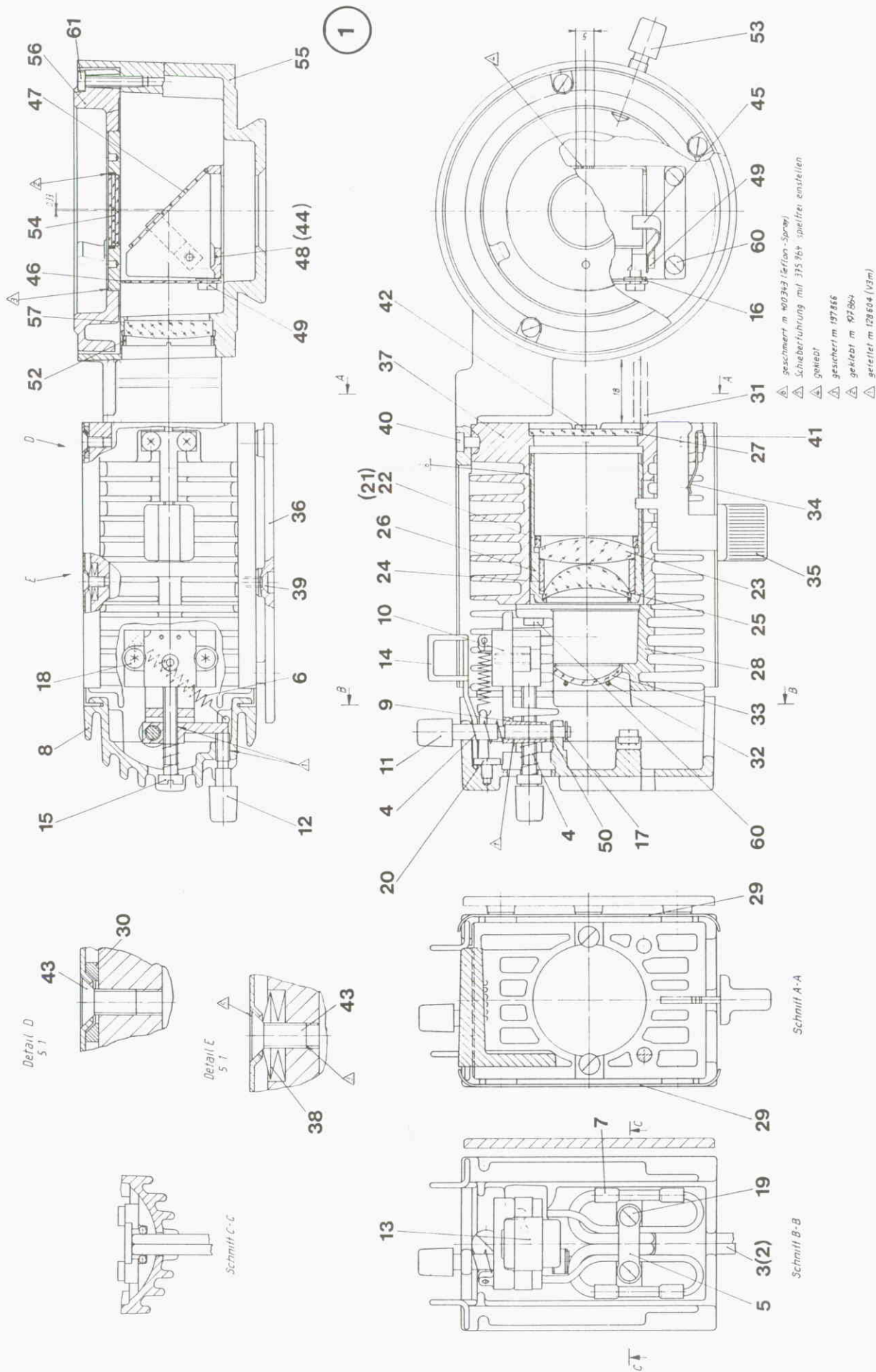
Attention:

The spares listet on this page have to be ordered directly from Leica AG, Heerbrugg with the 6-digit order No. of column " Bemerkung" (except for No 1 = order from Wetzlar).

Please address all inquiries regarding training, sevice documentation, warranty matters etc. for these products to Leica Heerbrugg AG.

Nr.	Bestell-Nr.	Stück	Benennung		Bemerkung
1	020-462.071-000	1	Auflichteinrichtung incident light attach.	dispositif-Epi dispositivo-Epi	V 599 007
2	- . .	1	Deckel verdrahtet cover wired	couvercle tapa	398 170
3	- . .	1	Kabel mit Stecker cable with plug	cable cable	160 821
4	- . .	1	Druckfeder pressure spring	ressort de traction muelle de traccion	250 216
5	- . .	1	Schelle clip	bride de serrage abrazadera	313 640
6	- . .	1	Zugfeder pull spring	ressort de traction muelle de traccion	336 615
7	- . .	1	Kabelstossverbinder cabel connector		343 783
8	- . .	1	Deckel cover	couvercle tapa	386 517
9	- . .	1	Lampenaufnahme lamp mount	monture de lampe montura para lampara	386 519
10	- . .	1	Lampenhalter lamp holder	support de lampe soporte p. lampara	386 521
11	- . .	1	Rändelschraube knurled screw	vis moletée tornillo moleteado	386 524
12	- . .	1	Rändelschraube knurled screw	vis moletée tornillo moleteado	386 527
13	- . .	1	Lampenfassung lamp mount	douille de lampe portalamparas	393 483
14	- . .	1	Verschlussbügel locking bracket		394 022
15	- . .	1	Schraube screw	vis tornillo	427 478
16	- . .	1	Scheibe Ø3,2 washer	rondelle arandela	177 975
17	- . .	1	Sicherungsscheibe Ø2,3 c-clip	disque de securite arandela de seguridad	243 211
18	- . .	1	Schraube screw	vis tornillo	309 763 M 3x10
19	- . .	1	Schraube screw	vis tornillo	348 985 M 3x8
20	- . .	1	Schraube screw	vis tornillo	358 316 M 3x4
21	- . .	1	Kollektor collector	collecteur colector	246 635
22	- . .	1	Gewinding threaded ring	bague filetee anillo con rosca	127 506
23	- . .	1	Kollektorlinse collector lens	collecteur colector	181 241
24	- . .	1	Distanzring spacer	bague d'entretoise anillo distanciador	394 023
25	- . .	1	Kollektorlinse collector lens	collecteur colector	398 881
26	- . .	1	Kollektorfassung collector mount		432 067
27	- . .	1	Wärmeschutzfilter heat absorption filter	filtre anti-calorique filtro anticálorico	126 119
28	- . .	1	Lampenkühler lamp cooler		386 561
29	- . .	1	Abdeckung cover	couverture cubierta	246 635
30	- . .	1	Isolerring insulating ring	bague isolante anillo aislante	386 651
31	- . .	1	Rändelstift knurled pin		389 139
32	- . .	1	Drahtbügel wire clamp		394 026

Nr.	Bestell-Nr.	Stück	Benennung		Bemerkung
33	- - -	1	Hohlspiegel reflector	miroir concave reflector	394 029
34	- - -	1	Feder spring	ressort muelle	398 843
35	- - -	1	Schieber slider	coulisseau corredera	412 002
36	- - -	1	Seitenplatte side cover		427 437
37	- - -	1	Lampengehäuse lamp housing		432 066
38	- - -	1	Tellerfeder saucer spring		169 912
39	- - -	1	Schraube screw	vis tornillo	251 684 M 3x10
40	- - -	1	Schraube screw	vis tornillo	296 953 M 3x8
41	- - -	1	Schraube screw	vis tornillo	372 446 M 3x5
42	- - -	1	Schraube screw	vis tornillo	398 981 M 2x4
43	- - -	1	Schraube screw	vis tornillo	375 964 M 3x6
44	- - -	1	Spiegelträger mirror carrier		398 364
45	- - -	1	Feder spring	ressort muelle	180 644
46	- - -	1	Polarisator polarizer	polarisateur polarizador	246 862
47	- - -	1	Strahlenteiler beam splitter		246 864
48	- - -	1	Spiegelträger mirror bracket	support de miroir portaespejos	400 125
49	- - -	1	Schraube screw	vis tornillo	175 868 M 3x6
50	- - -	1	Scheibe Ø3,2 washer	rondelle arandela	177 975
51	- - -	1	Niedervoltlampe low voltage lamp		398 365
52	- - -	1	Gewinding threaded ring	bage filetee anillo con rosca	127 506
53	- - -	1	Klemmschraube clamping screw		377 516
54	- - -	1	Zirkularpolfilter circular pol. filter		398 363
55	- - -	1	Gehäuse housing	boitier caja	400 122
56	- - -	1	Tubusträger tube carrier		400 124
57	- - -	1	Beleuchtungslinse illuminating lens	lentille d'eclairage lente de iluminacion	410 211
58	- - -	1	Schutzdeckel protective cap		424 316
59	- - -	1	Schutzdeckel protective cap		327 436
60	- - -	1	Schraube screw	vis tornillo	296 953 M 3x8
61	- - -	1	Schraube screw	vis tornillo	381 722 M 3x16

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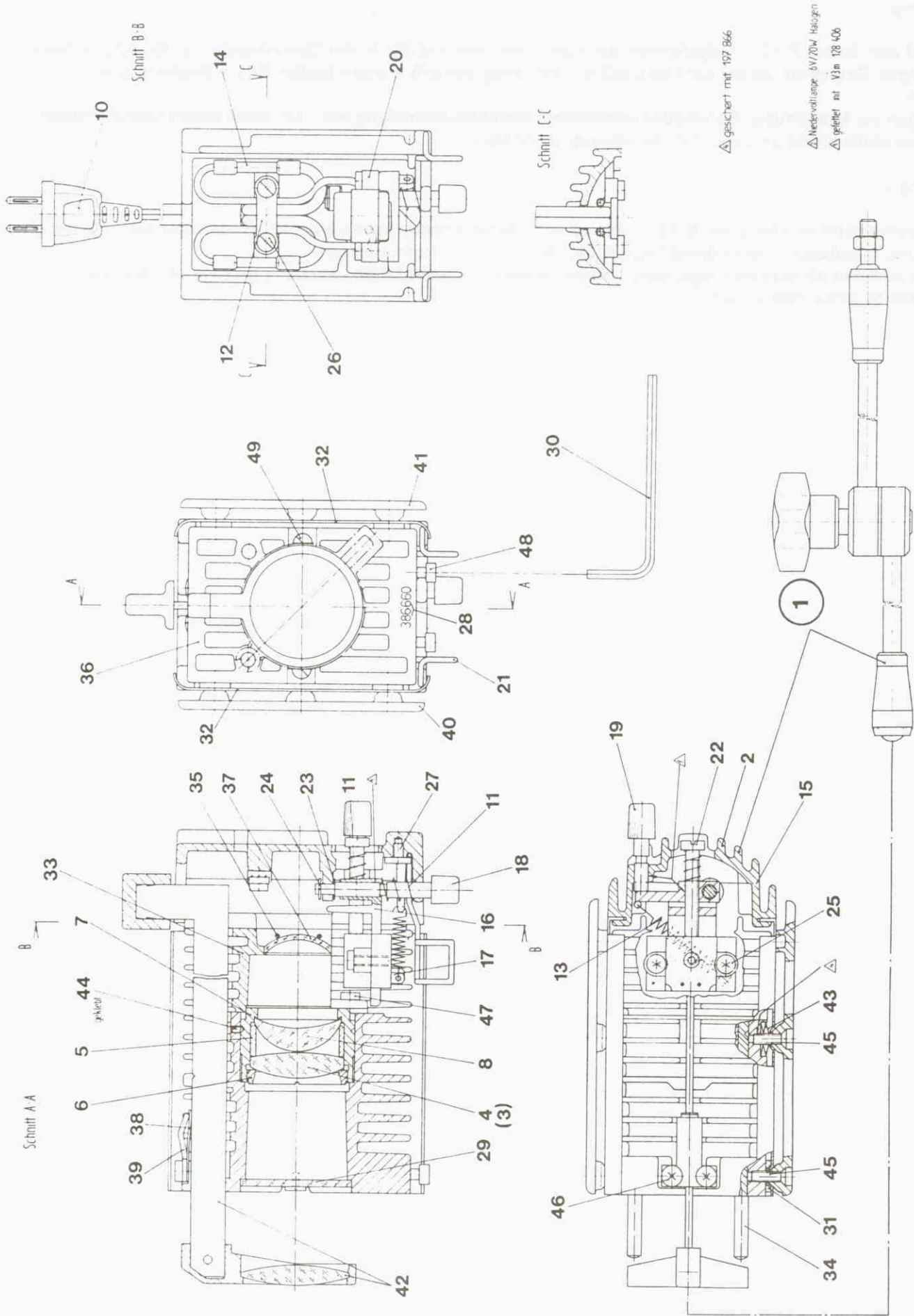
Attention:

The spares listet on this page B 12.1 have to be ordered directly from Leica AG, Heerbrugg with the 6-digit order No. in column " Bemerkung" (except for No 1 = order from Wetzlar).

Please address all inquiries regarding training, service documentation, warranty matters etc. for these products to Leica Heerbrugg AG.

Nr.	Bestell-Nr.	Stück	Benennung		Bemerkung
1	020-462.073-000	1	Lampenhaus mit Gelenkarm lamp housing	boitier de lampe caja de lampara	V 599 009
2	020-462.016-021	1	Lampenhaus lamp housing	boitier de lampe caja de lampara	
3	- . .	1	Kollektor collector	collecteur colector	398 169
4	- . .	1	Kollektorlinse collector lens		181 241
5	- . .	1	Distanzring spacer	bague d'entretoise anillo distanciador	394 023
6	- . .	1	Gewinding threaded ring	bague filetee anillo con rosca	394 025
7	- . .	1	Kollektorfassung collector mount		398 880
8	- . .	1	Kollektorlinse collector lens		398 881
9	- . .	1	Deckel cover	couvercle tapa	398 170
10	- . .	1	Kabel mit Stecker cable with plug	cable cable	160 821
11	- . .	1	Druckfeder pressure spring	ressort de pression muelle de compresion	250 216
12	- . .	1	Schelle clip	bride de serrage abrazadera	313 640
13	- . .	1	Zugfeder pull spring	ressort de traction muelle de traccion	336 615
14	- . .	1	Kabelstossverbinder cable connector		343 783
15	- . .	1	Deckel cover	couvercle tapa	386 517
16	- . .	1	Lampenführung lamp guide		386 519
17	- . .	1	Lampenträger lamp carrier		386 521
18	- . .	1	Rändelschraube knurled screw	vis moletée tornillo moleteado	386 524
19	- . .	1	Rändelschraube knurled screw	vis moletée tornillo moleteado	386 527
20	- . .	1	Lampenfassung lamp mount	douille de lampe portalamparas	393 483
21	- . .	1	Verschlussbügel locking bracket		394 022
22	- . .	1	Schraube screw	vis tornillo	427 478
23	- . .	1	Scheibe Ø 3,2 washer	rondelle arandela	177 975
24	- . .	1	Sicherungsscheibe c-clip	disque de securite arandela de seguridad	243 211
25	- . .	1	Schraube screw	vis tornillo	309 763 M 3x10
26	- . .	1	Schraube screw	vis tornillo	348 985 M 3x8
27	- . .	1	Schraube screw	vis tornillo	358 316 M 3x4
28	- . .	1	Niedervoltlampe low voltage lamp		386 660
29	- . .	1	Wärmeschutzfilter heat absorption filter	filtre anti-calorique filtro anticalorico	126 119
30	- . .	1	Inbusstiftschlüssel allen key		229 341
31	- . .	1	Lampenkühler lamp cooler		386 561

Nr.	Bestell-Nr.	Stück	Benennung		Bemerkung
32	- - -	1	Abdeckung cover	couverture cubierta	386 562
33	- - -	1	Isolerring insulating ring	bague isolante anillo aislante	386 651
34	- - -	1	Rändelstift knurled pin		389 139
35	- - -	1	Drahtbügel wire clamp		394 026
36	- - -	1	Lampengehäuse lamp housing		394 028
37	- - -	1	Hohlspiegel reflector	miroir concave reflector	394 029
38	- - -	1	Feder spring	ressort muelle	398 843
39	- - -	1	Bride cable clip		421 828
40	- - -	1	Seitenplatte side plate		427 437
41	- - -	1	Seitenplatte beschriftet side plate		427 438
42	- - -	1	Schieber mit Beleuchtungslinse slide with illumination lens		438 111
43	- - -	1	Tellerfeder saucer spring		169 912
44	- - -	1	Gewindestift threaded pin	goupille filetee perno con rosca	176 091 M 2x3
45	- - -	1	Schraube screw	vis tornillo	251 684 M 3x10
46	- - -	1	Schraube screw	vis tornillo	276 159 M 3x6
47	- - -	1	Schraube screw	vis tornillo	296 953 M 3x8
48	- - -	1	Schraube screw	vis tornillo	309 763 M 3x10
49	- - -	1	Schraube screw	vis tornillo	398 981 M 2x4

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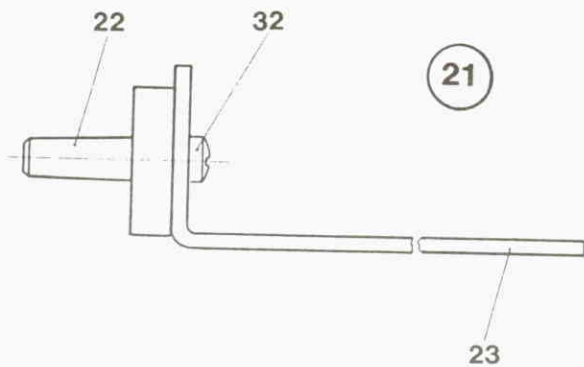
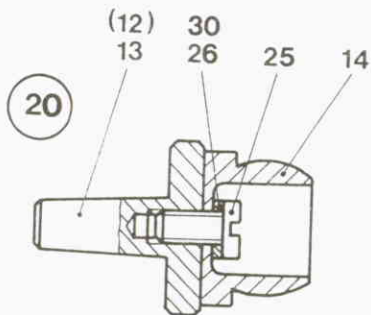
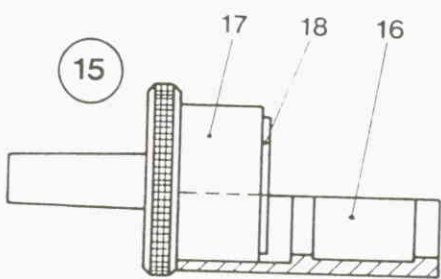
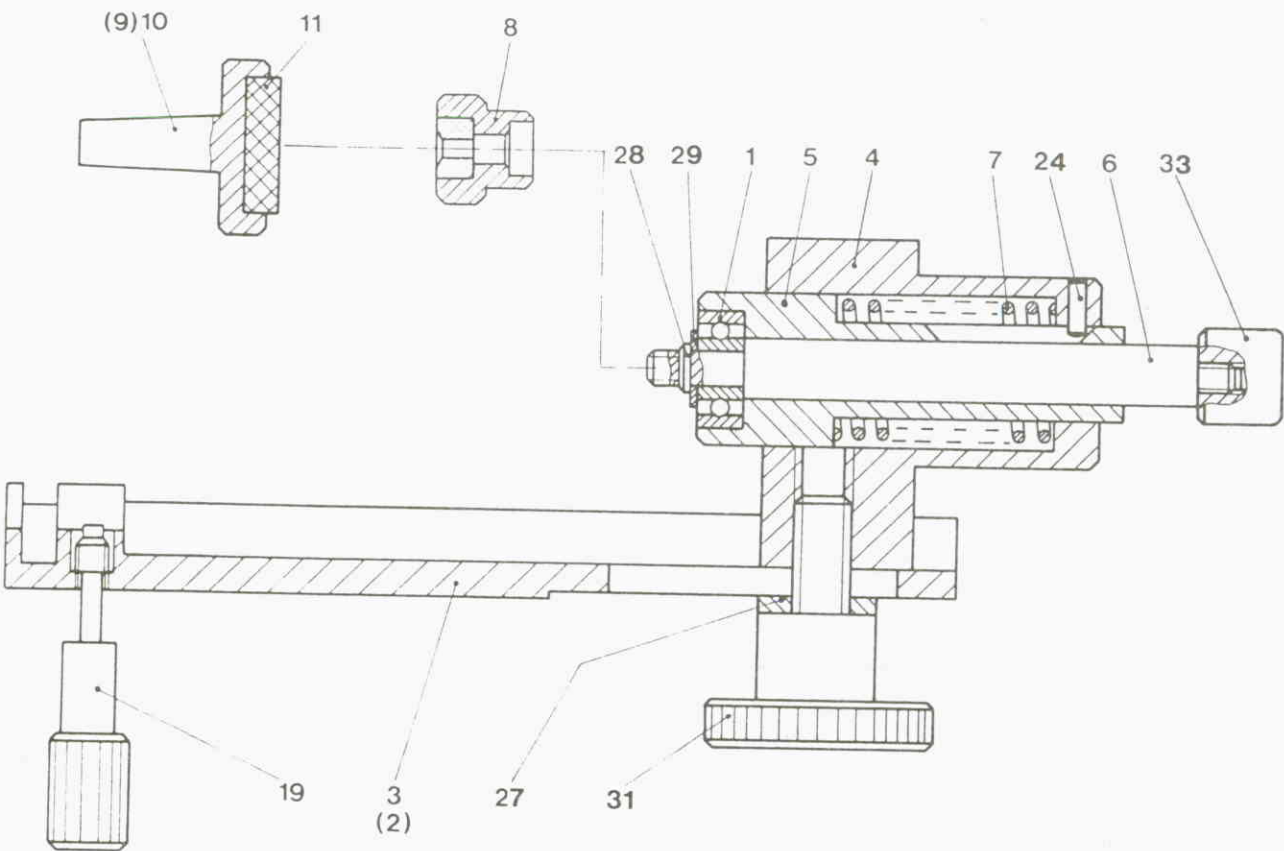
Attention:








The spares listet on page B 13.1 have to be ordered directly from Leica AG, Heerbrugg with the 6-digit order No. of column " Bemerkung" (except for No 1 and 2 = order from Wetzlar).

Please address all inquiries regarding training, sevice documentation, warranty matters etc. for these products to Leica Heerbrugg AG.

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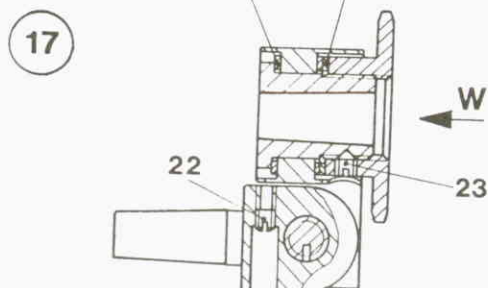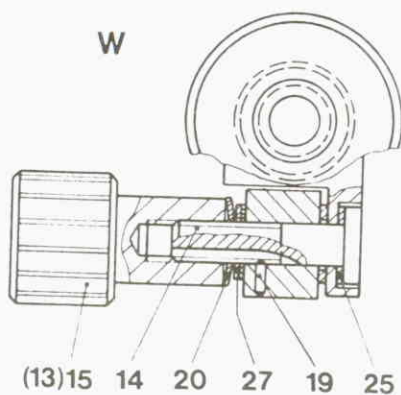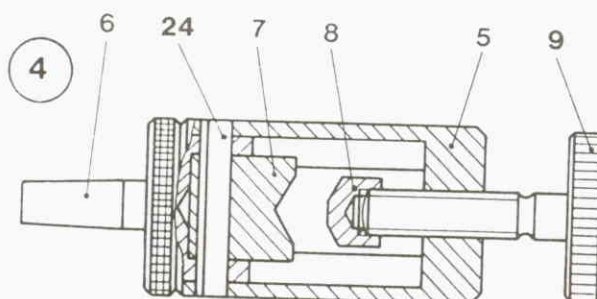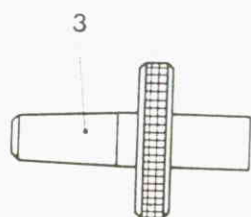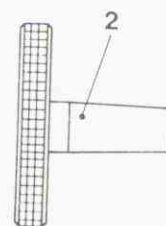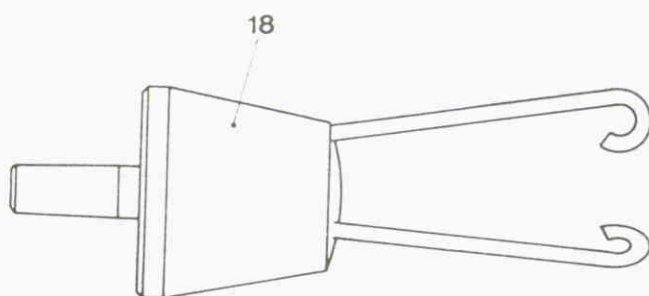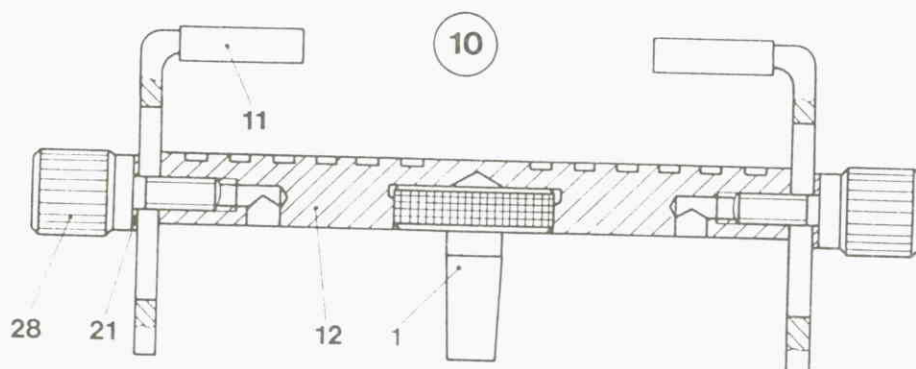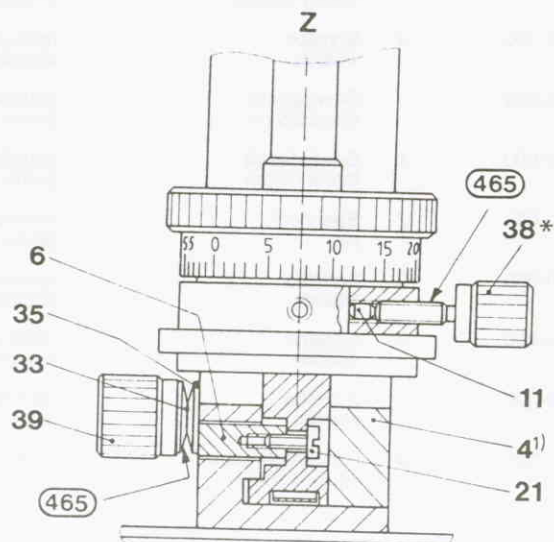
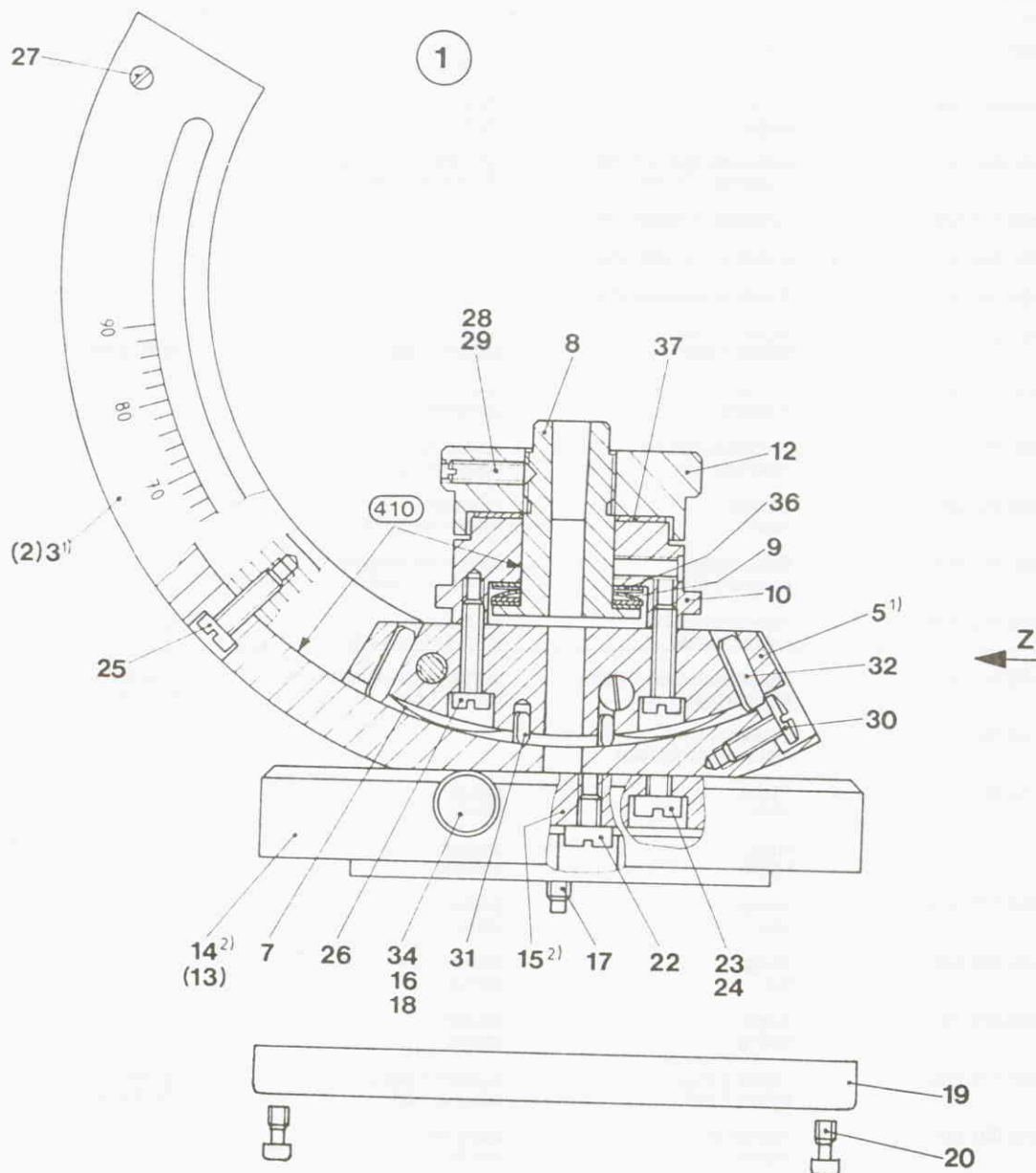
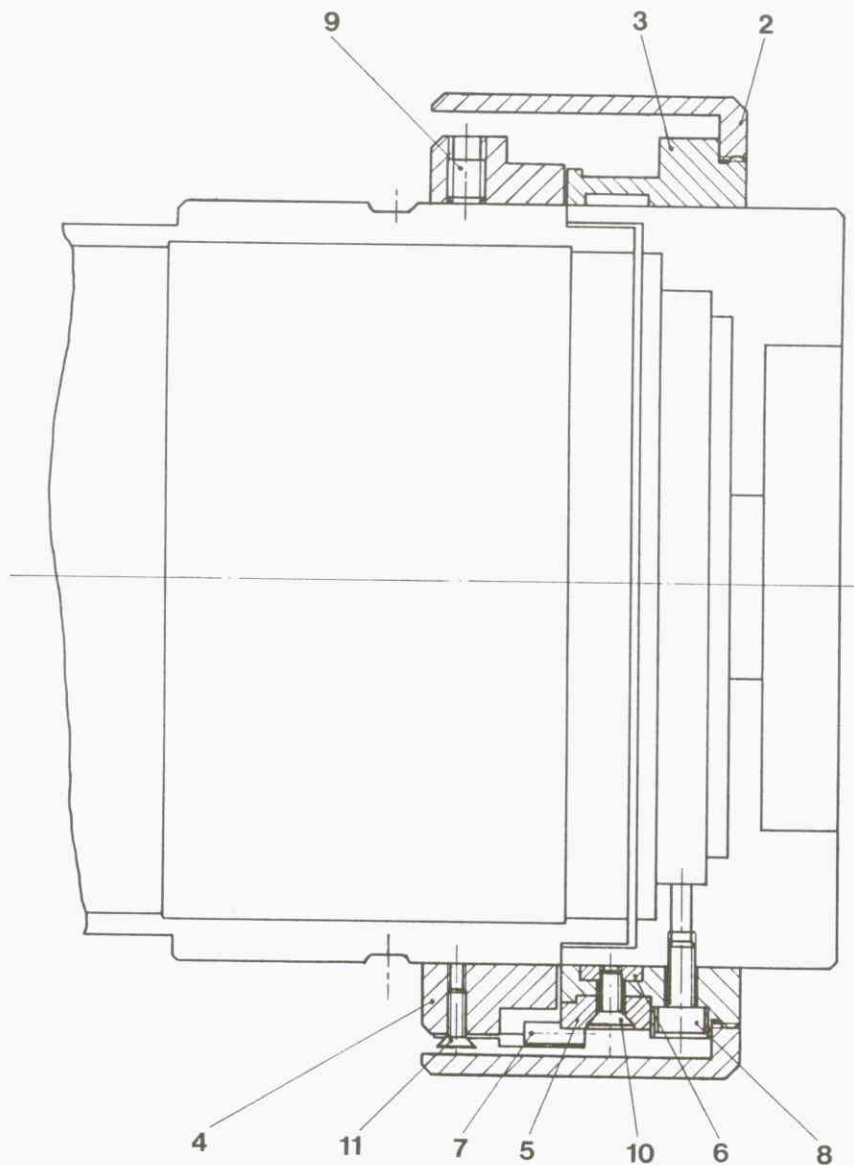
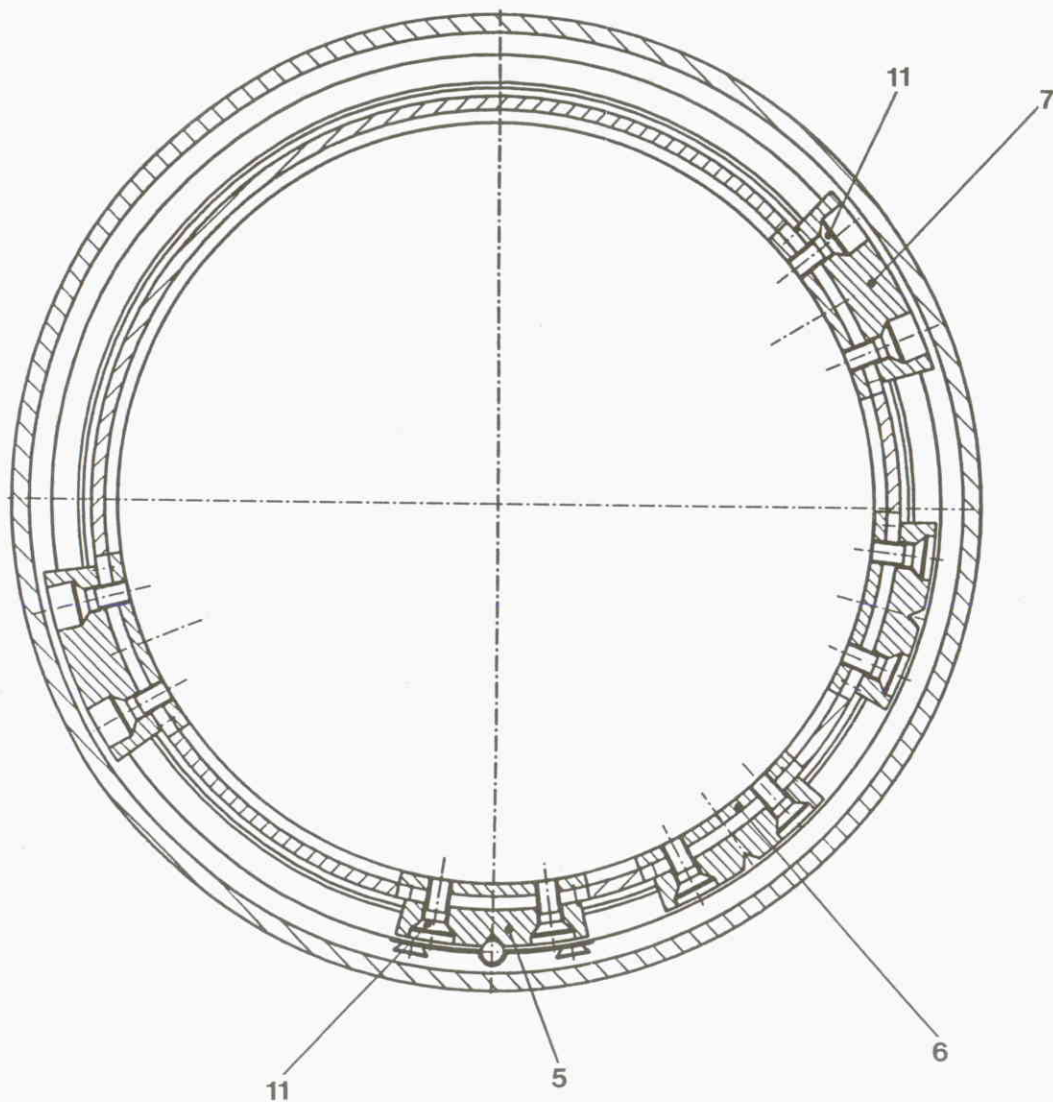
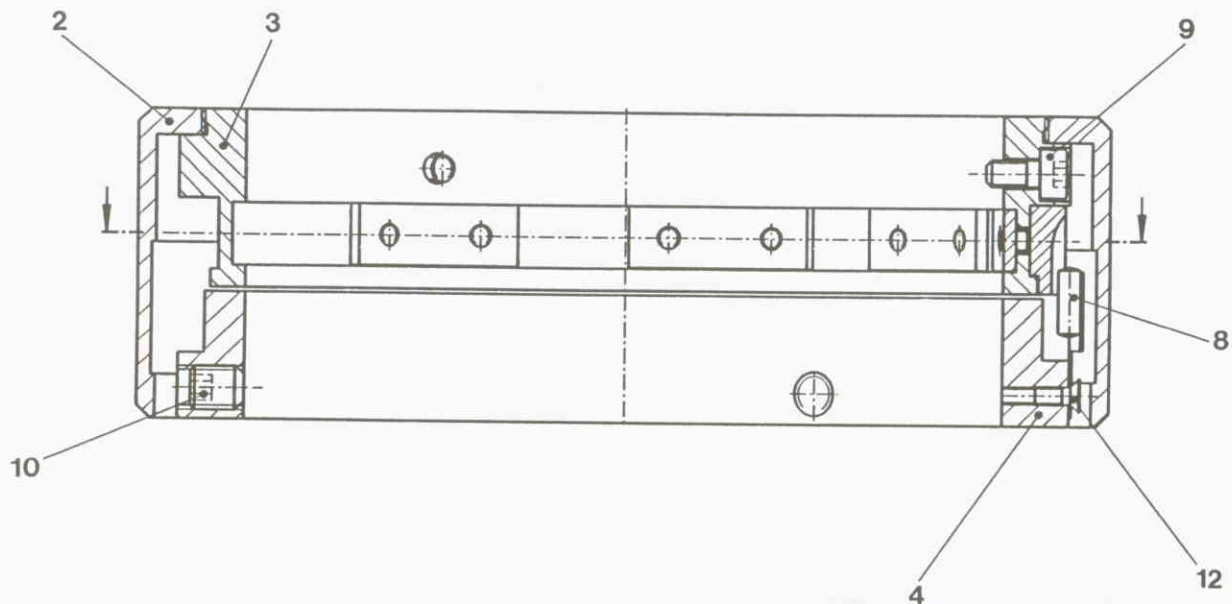
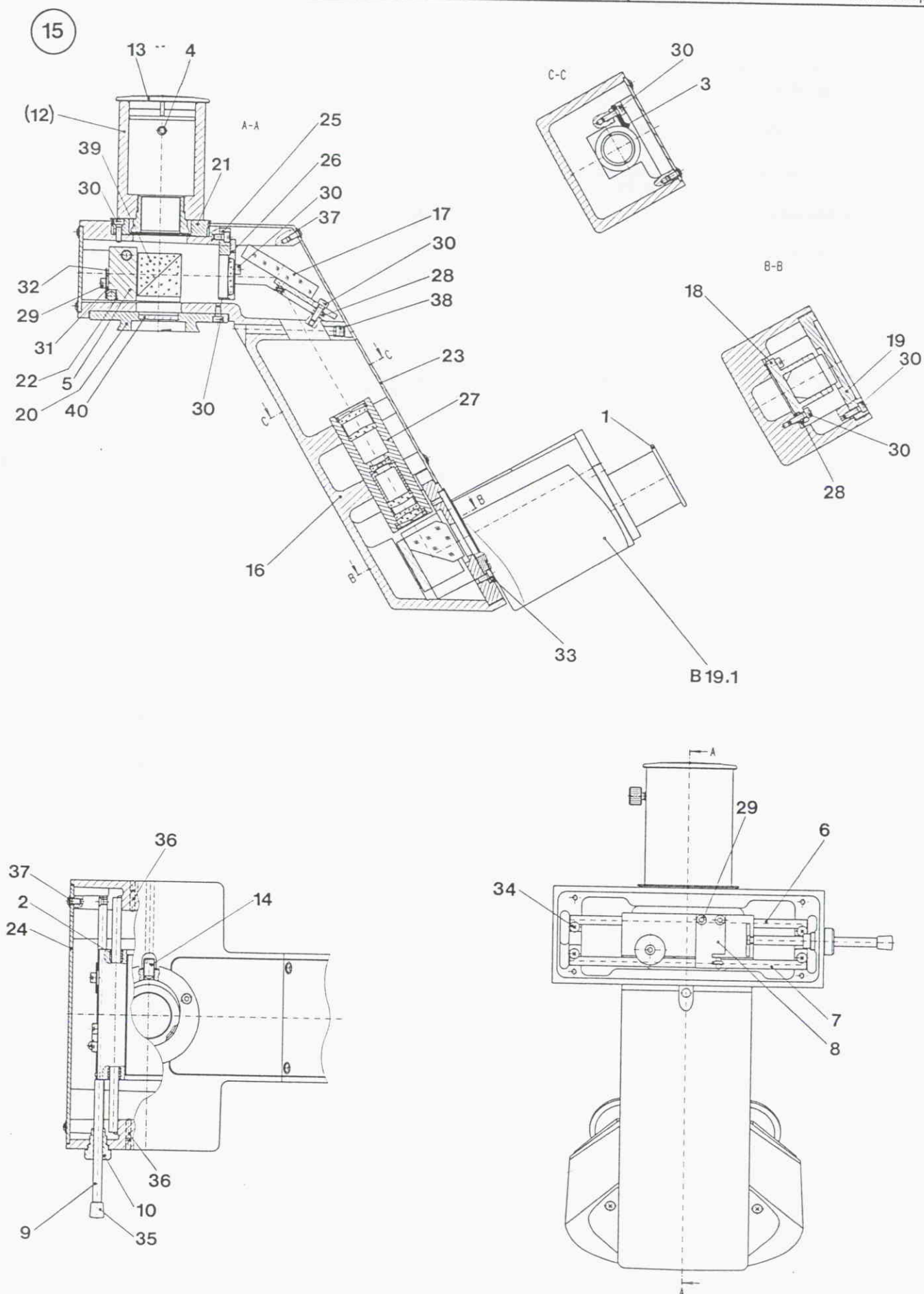
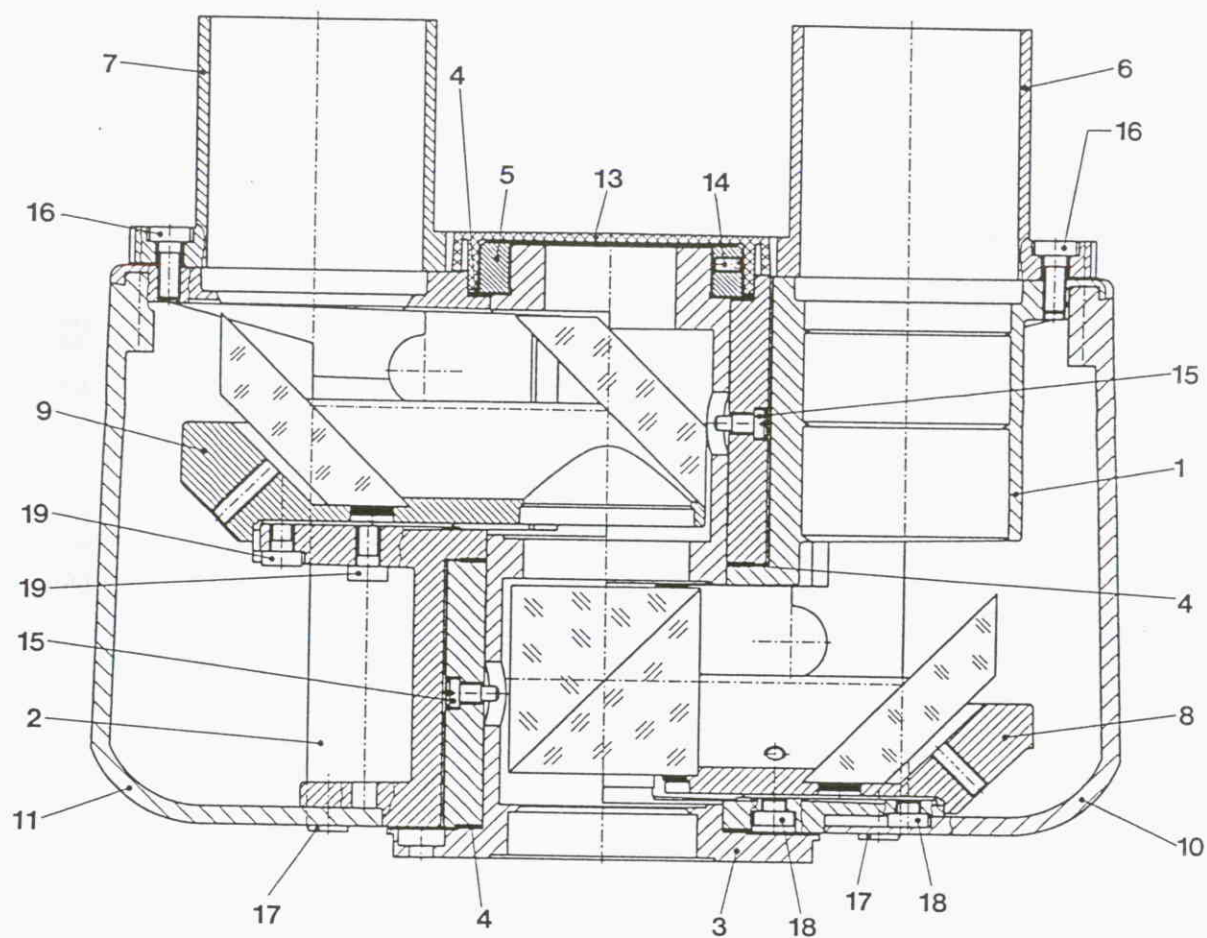
Disassembly of Z- drive, right-hand side

Unscrew the 4 screws (2) of the lid (1).
Slacken the screw (3) by about 3-4 rotations. Carefully pull out the axle (4), then remove the washer (5), gearwheel (6) and sleeve (7) one after the other.
Lay the microscope on its back and unscrew the base plate. Move the x adjustment to the centre position so that the screw holes of the coarse and fine drive coincide. Then unscrew the 2 screws (8) of the coarse and fine drive. Pull the drive out towards the right.
Slacken the screw (9) by about 3-4 rotations. Carefully pull out the axle (10), then remove the washer (11) and sleeve (12) one after the other. Then unscrew the 2 screws (14) holding the holder with the gearwheel (13). Slacken the 4 pressure screws (15) by about 5 rotations.
Unscrew the 4 screws (16). Remove the strip (17) with the stage bracket (18), roller cages (19) and rollers (20).
Clean all parts and grease with the specified lubricant. The components are reassembled in the reverse order.

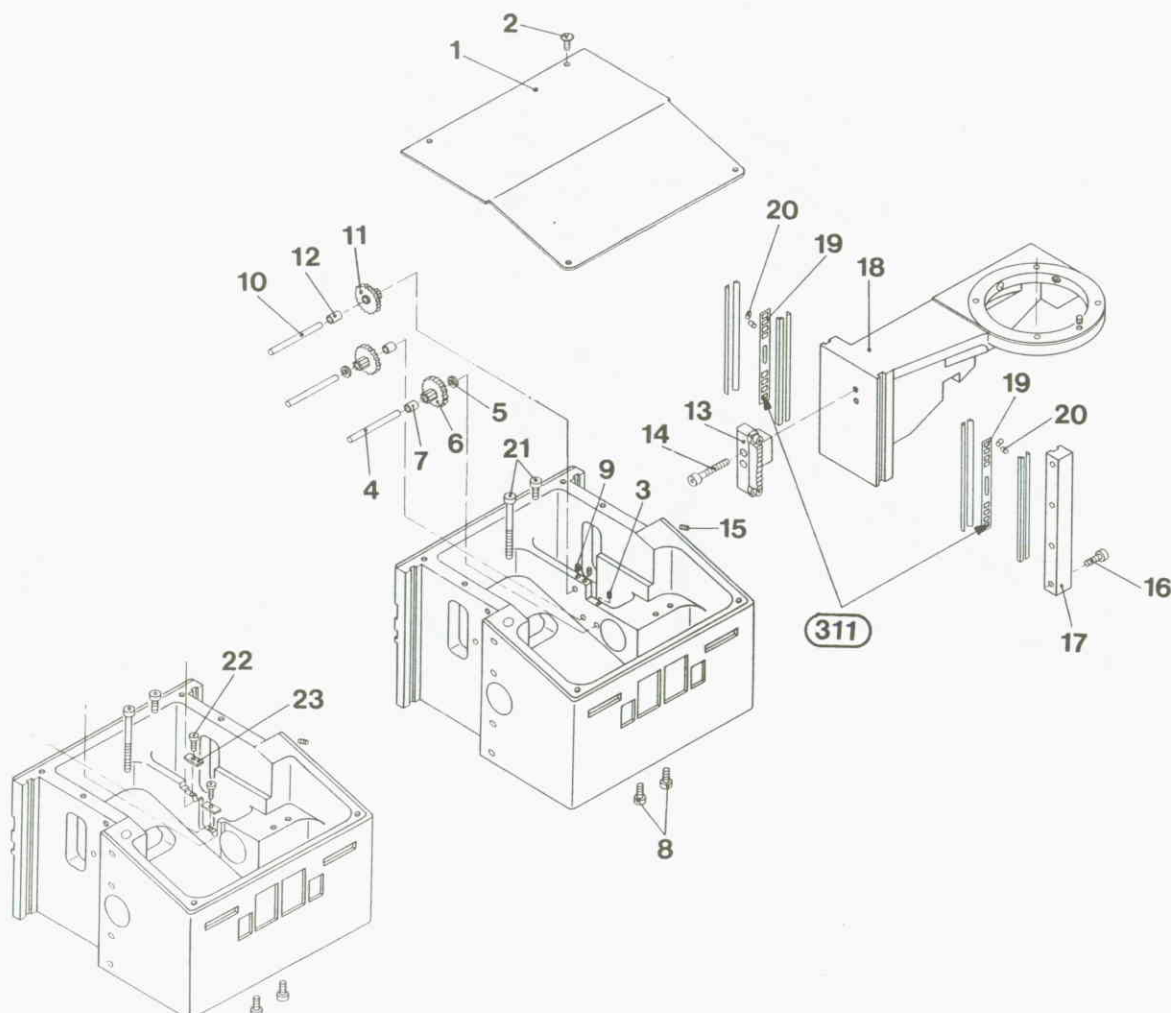
Disassembly of Z drive, left-hand side

The disassembly of the Z drive on the left-hand side is done in the same order, except that for the removal of the coarse and fine adjustment the 2 screws (21) have to be removed instead of the screws (8), and the microscope need not be laid on its back.

Attention !

The mounting of the axles (10) with the gearwheels (11) was modified from serial number 176897, construction date 11.96.

Before the gearwheels (11) can be disassembled, the screws (22) have to be unscrewed and then the retaining discs (23) removed.

Disassembly of x -adjustment

It is easier to disassemble or assemble the x adjustment if the two z adjustments on the left- and right-hand side are disassembled first. See page D 1.1.

Screw off the lid (1) with 4 screws (2).

Make a sketch of the wiring and write down the colours of the wires.

Pull the wires off the switches (3).

Unscrew the clamp fitting (4) with the screw (5). Unscrew the 2 screws (6) with the holders (7) and potentiometer (11) for light adjustment, then slacken the screw (8) and remove the graduated disk (9). Screw off the nut (10) and take the potentiometer (11) out of its holder (7).

Slacken the screw (12) in the knurled knob (13) of the x adjustment. Pull off the knob (13). Unscrew the 2 locknuts (14).

Unscrew the two screws (15) and take the spindle (16) with spindle nut (17) (upend) and the washer (27) out of the casing towards the left.

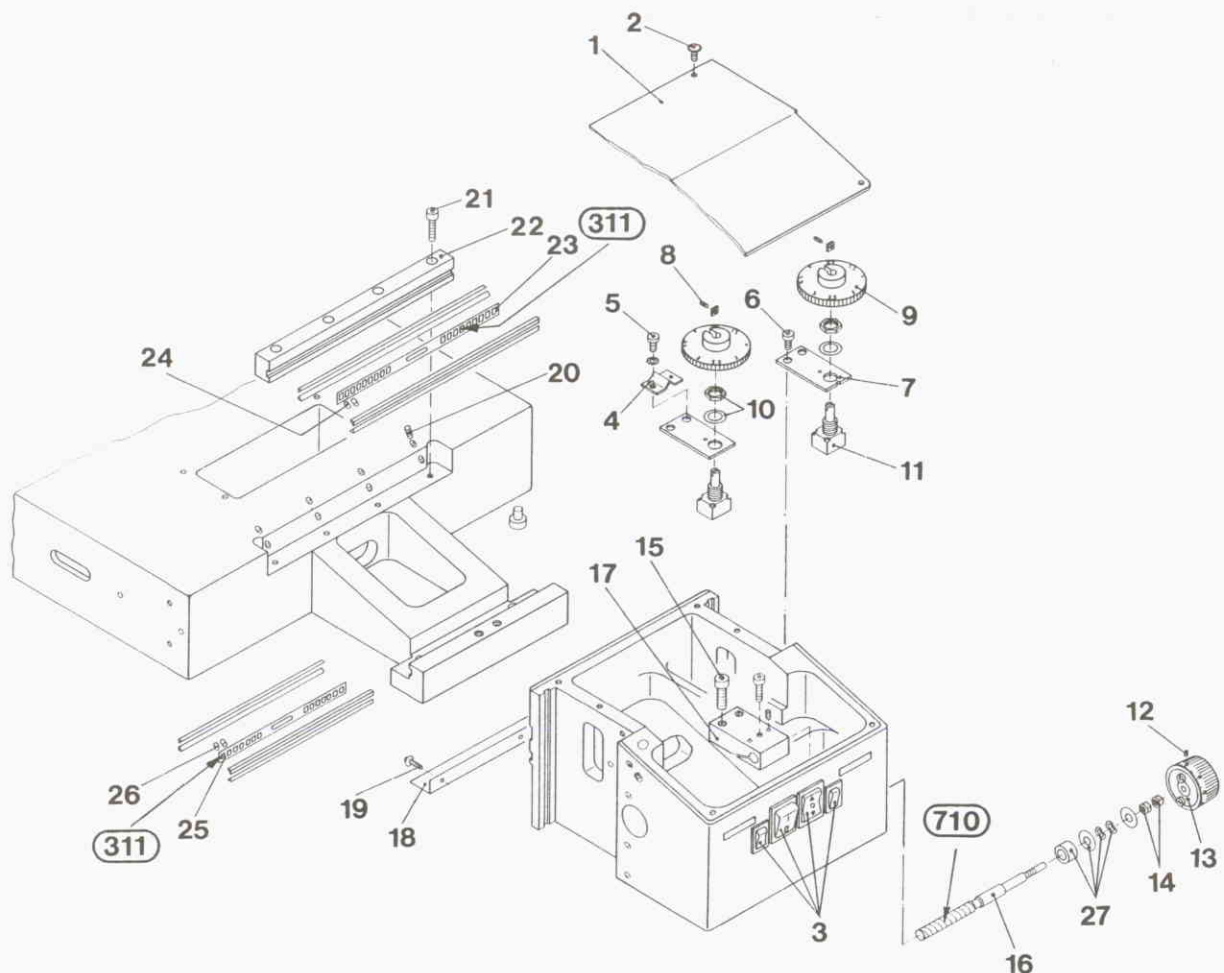
Unscrew the cover (18) with 3 screws (19).

Slacken 4 pressure screws (20) by about 5 rotations.

Remove 4 screws (21) and take out strip (22) with roller cage (23) and rollers (24).


Push the gearbox to the back, pick up from the front and lift off upwards. Be careful not to drop the roller cage (25) and the rollers (26). Clean all parts and grease with specified lubricants.

The components are reassembled in the reverse

Disassembly of column with motor (z height adjustment)




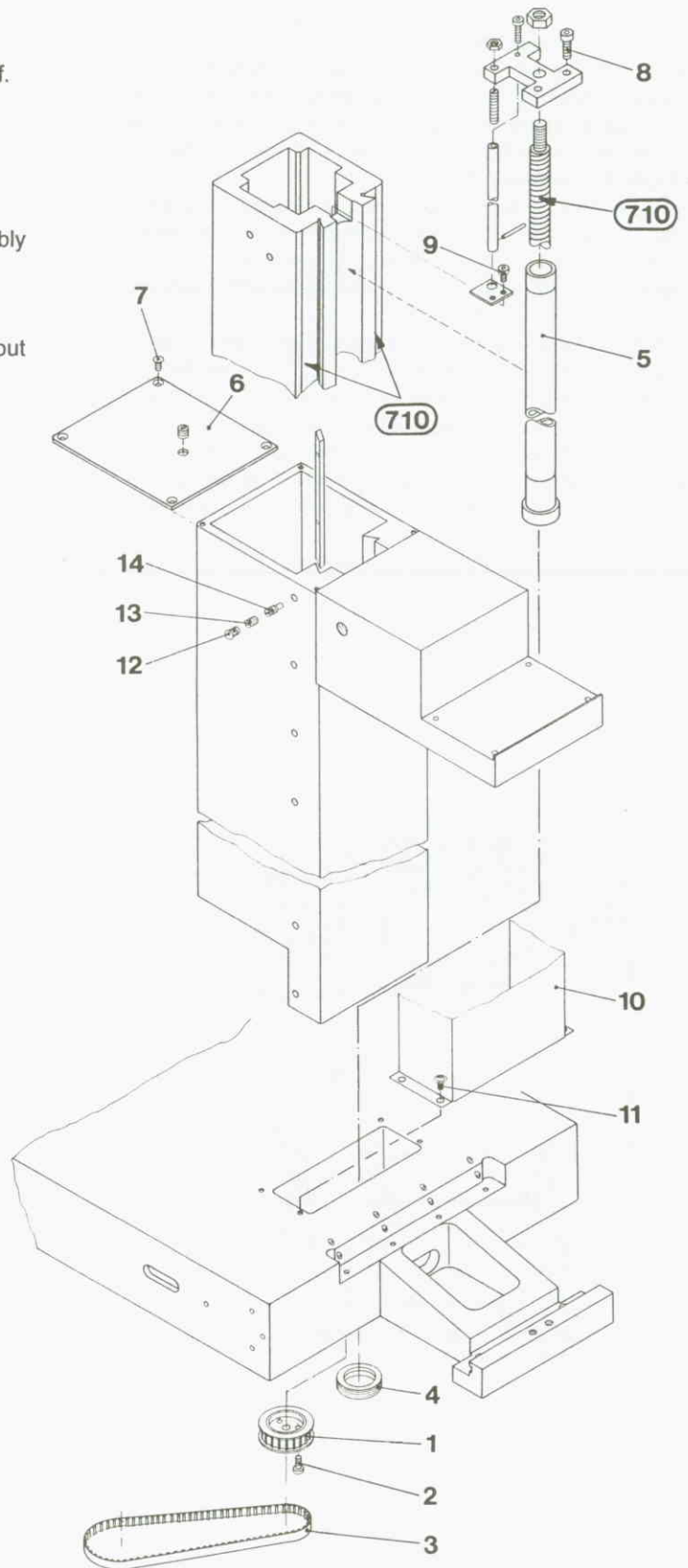
Screw off the comparison bridge with 4 screws.
Lay the microscope on its back and screw off the base plate.
Unscrew the crown gear (1) with 3 screws (2) and lift off.
Remove the toothed belt (3).
Unscrew the ring nut (4) and unscrew the tube (5) downwards.
Screw off the lid (6) with 4 screws (7).
Unscrew 2 screws (8).
Unscrew 2 screws (9), then lift out the whole subassembly with the spiral spindle.
Screw off the cover plate (10) with 4 screws (11).
Remove 7 caps (12), then unscrew 7 screws (13) and slacken 7 screws (14) by about 2 rotations. Then push out the outer column with support upwards.
Clean all parts and grease with the specified lubricants.
The components are reassembled in the reverse order.

Assembly of zoom device

Screw the two screws (1) of the zoom objective in as far as the stop and then screw back by $\frac{1}{2}$ rotation. Push the ring (2) over the objective from above (catch spring points upwards). Then push the clickstop ring (3) (clickstops point downwards) over the objective from above until the two drill holes of screw (4) coincide with the holes of screw (1). The ring (3) is secured in the screw holes of (1) with the two M 3x6 DIN 912 screws (4).

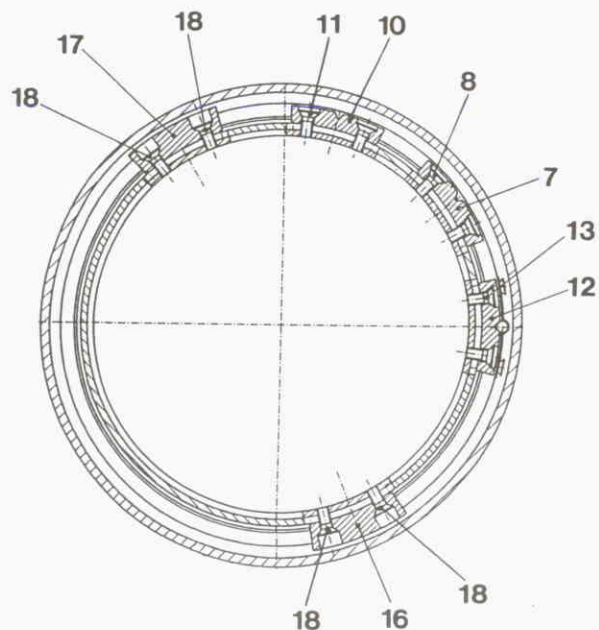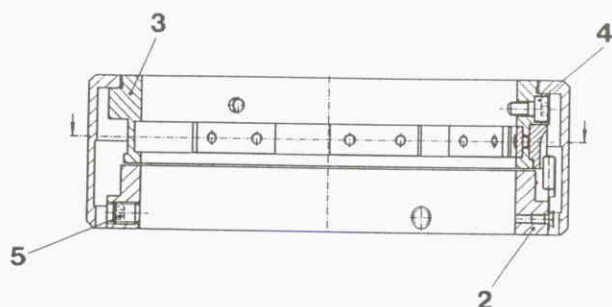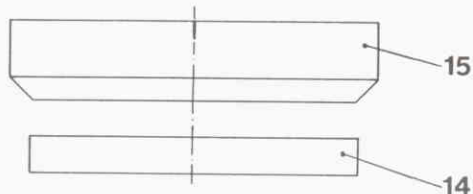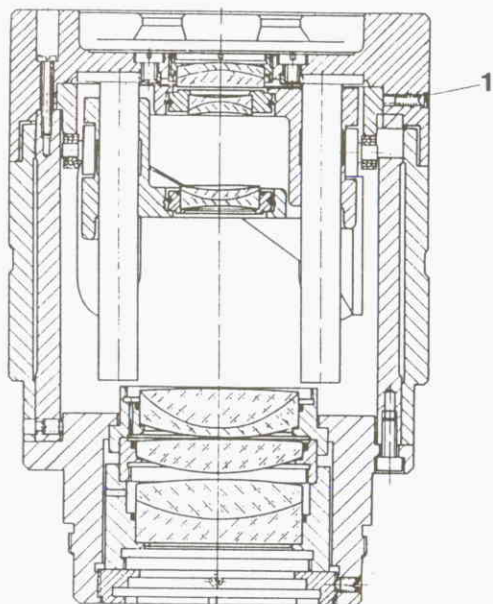On the left-hand objective, push the lower ring (2) against the ring (3). Set the Makrozoom objective to magnification 12.5x, then engage the catch spring (6) and the middle clickstop (7) in this position and screw down with screws (8 and 9). Then tighten the lower ring (2) slightly with three M 5x6 DIN 913 screws (5). The distance between the two rings (2 and 3) should be between 0.1 and 0.2 mm. Now tighten the screws by an equal amount, making sure that the objective can still be easily rotated. Now set the objective magnification to 8x, engage the right-hand clickstop (10) under the catch spring (6) and screw tight with the screws (11). Now set the objective magnification to 20x, engage the left-hand clickstop (12) under the catch spring (6) and screw tight with the screws (13).

Adjustment: Right-hand objective

Screw the quarter-wave compensator off the objective, if applicable. Screw the ring (14) off the objective, then push a new index ring (15) onto the objective from the bottom and tighten slightly with the ring (14). Mount the objectives onto the comparison bridge and screw on with screw (16). Put stage micrometers on both stages and form an image with the split image setting of the comparison bridge. Align the scale marks of the stage micrometers along the dividing line between the two part images as shown in Fig. 1. Set the right-hand objective at magnification 12.5x. Engage the catch spring (6) in clickstop (12). Adjust the clickstop (12) by sliding to the right and left until the two stage micrometers are imaged in the same scale ratio. Then fix the clickstop position (12) with the screws (13). The clickstop positions for 8x and 20x are set in the same way.

Adjustment: Upper and lower stop

The upper and lower stops for the two objectives are determined by sliding the stops (16 and 17) towards each other until the two stage micrometers (Fig. 1) are imaged in the same scale ratio.

The stops (16 and 17) are then fixed in position with the screws (18).

Set the objective to magnification 12.5x. Make the graduation mark of the index ring (15) coincide with the 12.5x mark of the Makrozoom objective, the ring (14) is then turned firmly against it. Lightly lubricate the clickstops with grease 410/01. Unscrew both objectives from the comparison bridge with screw (16). Screw the cover (19) over the clickstop device from the top. Screw the quarter-wave compensator back onto the right-hand objective if appropriate. Mount the objectives back onto the comparison bridge and secure with screw (16). Check the adjustment.

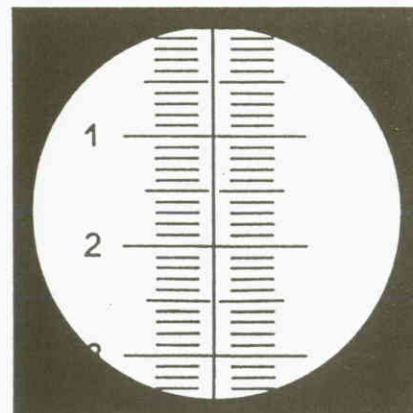
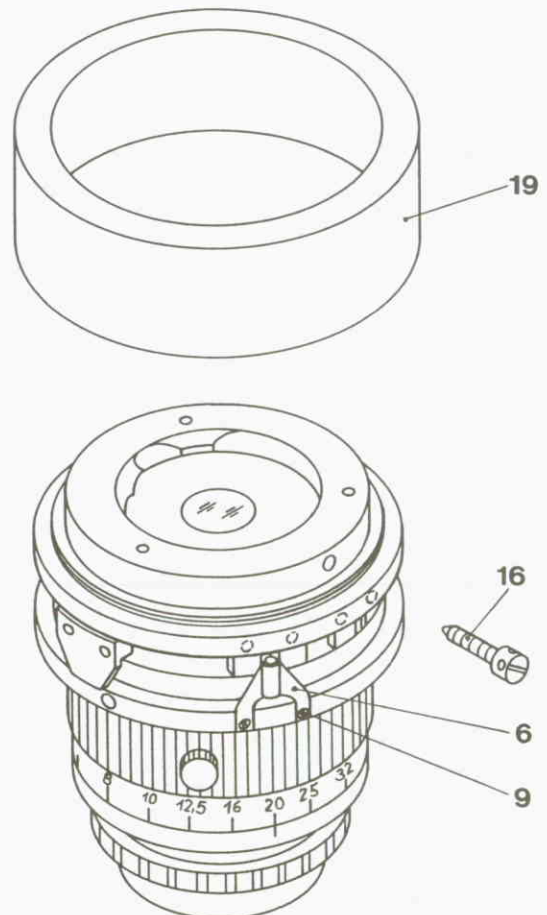
Fig. 1

The construction status of the microscope is indicated with a status number. This can be seen on the type plate on the back of the microscope after the letters BZ.

Construction status 00

Index number :

165 750 to 165 755

Construction status 01

Index number :

170 023 to 170 026
170 028 to 170 035
170 037 to 170 039

New additional transformer for the two 45 VA lamps.
See sheet B 4.2, item 2

Reason: elimination of brightness fall-off while
operating the Z height adjustment of the microscope.

Construction status 02

Index number:

171 466 to 171 473
171 477 to 171 506
176 816 to 176 908
193 577 to 193 533

Analog control 2x30 W power supply. See sheet B 4.2, item 6

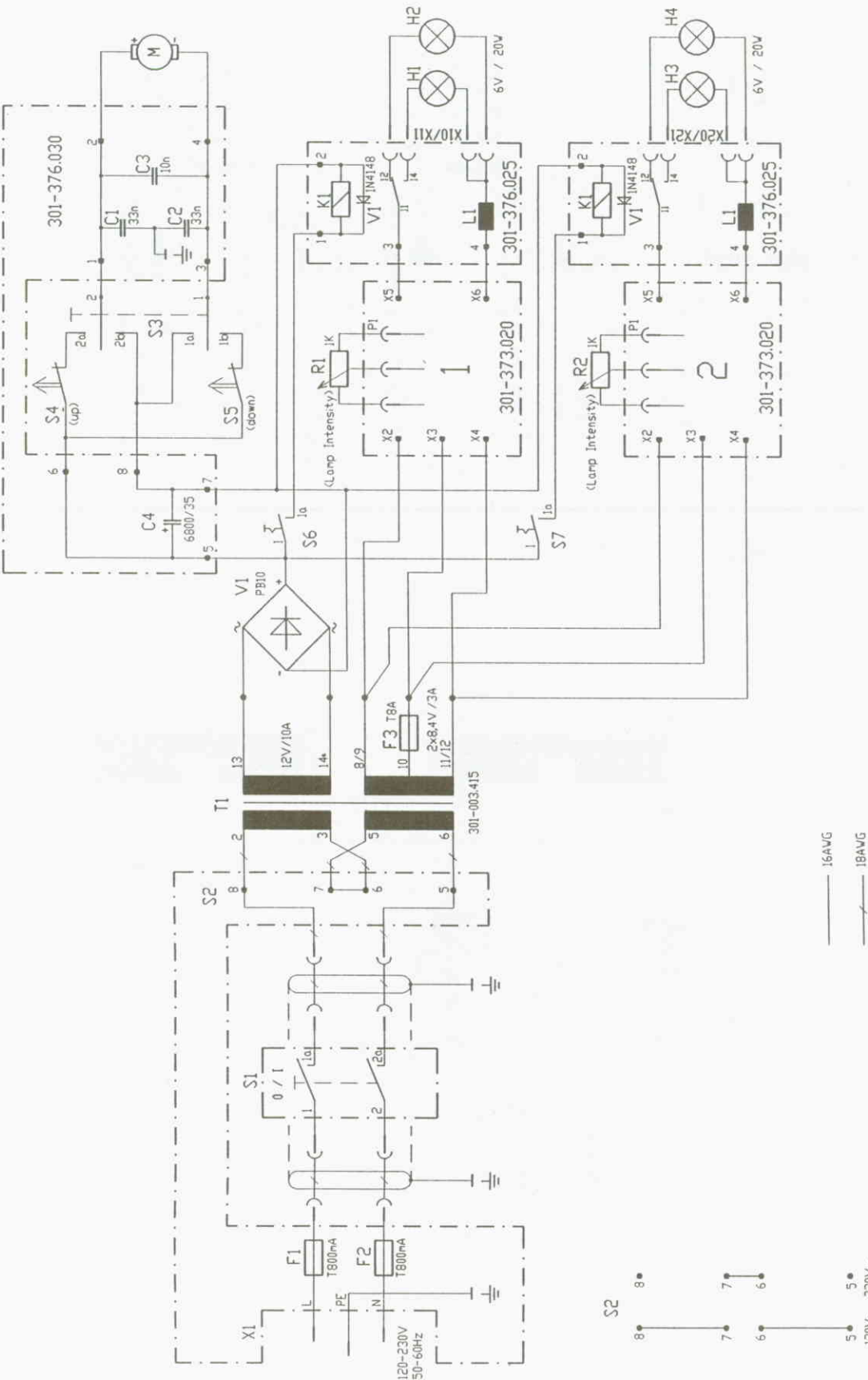
Reason: standardization of DM C board with that of DM IL

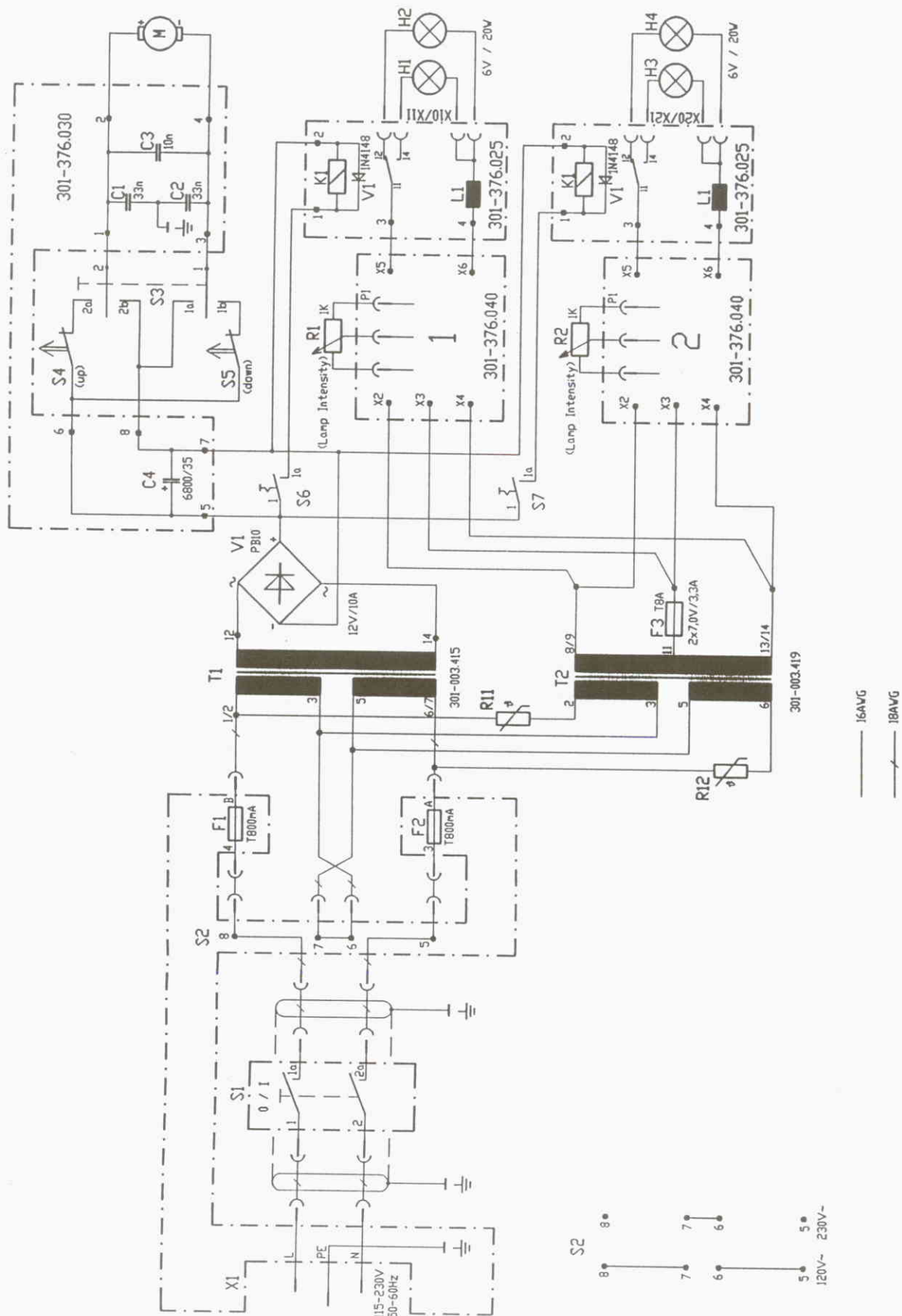
Construction status 03

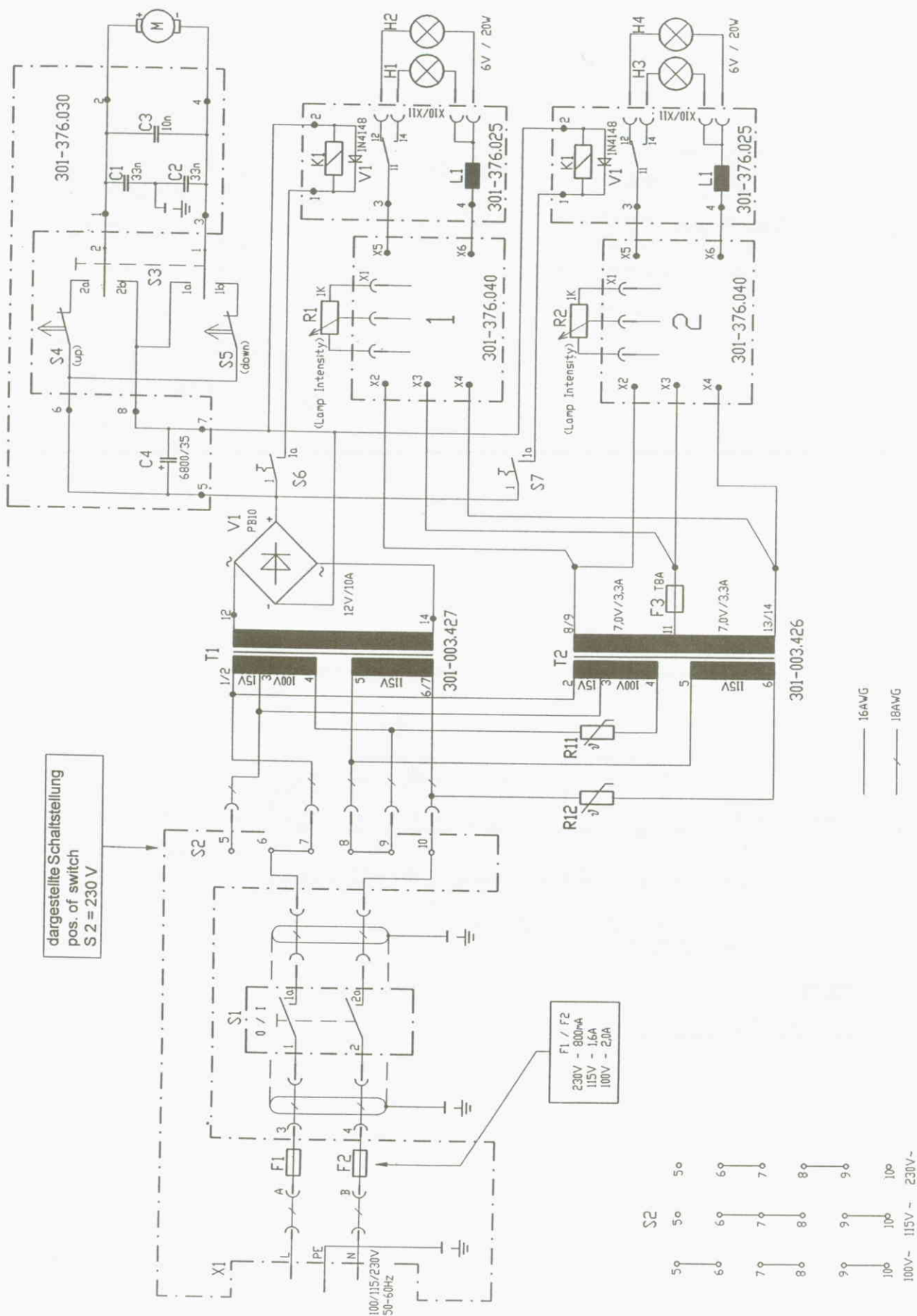
Index number:

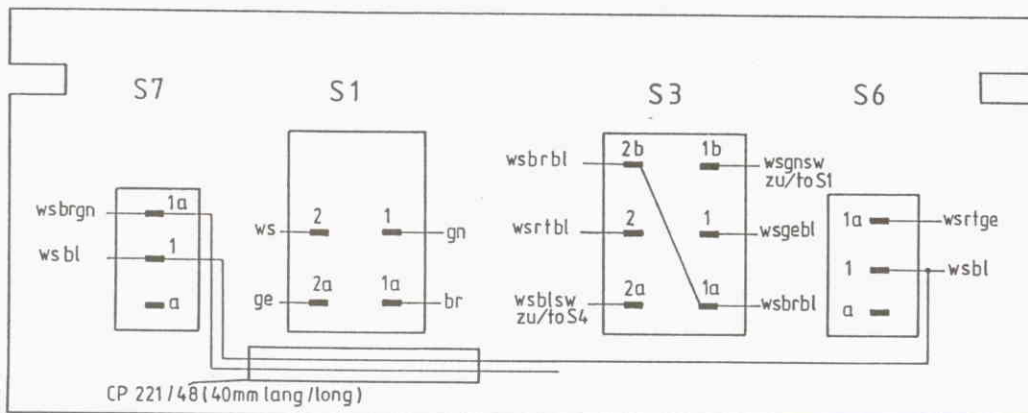
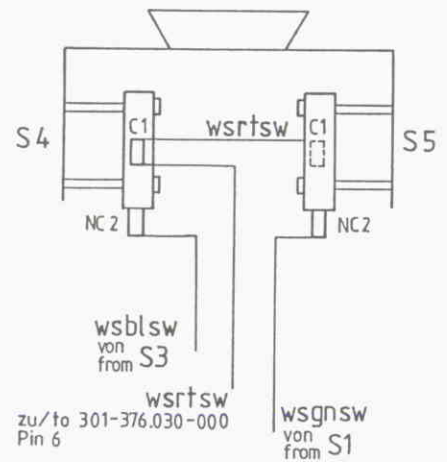
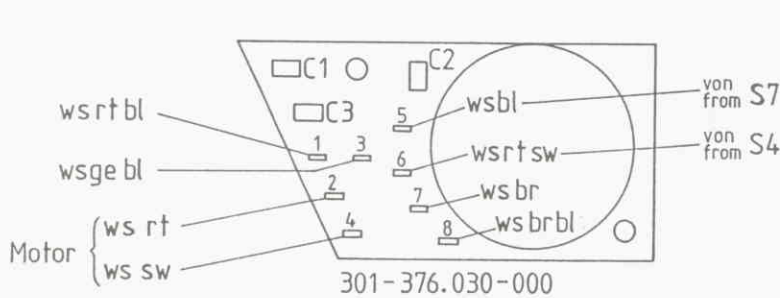
195 710 to --->

Use of transistors of only one manufacturer.
Reason: difference in light intensity between right- and left-hand
side.
See Service News MQM 03 FY 98

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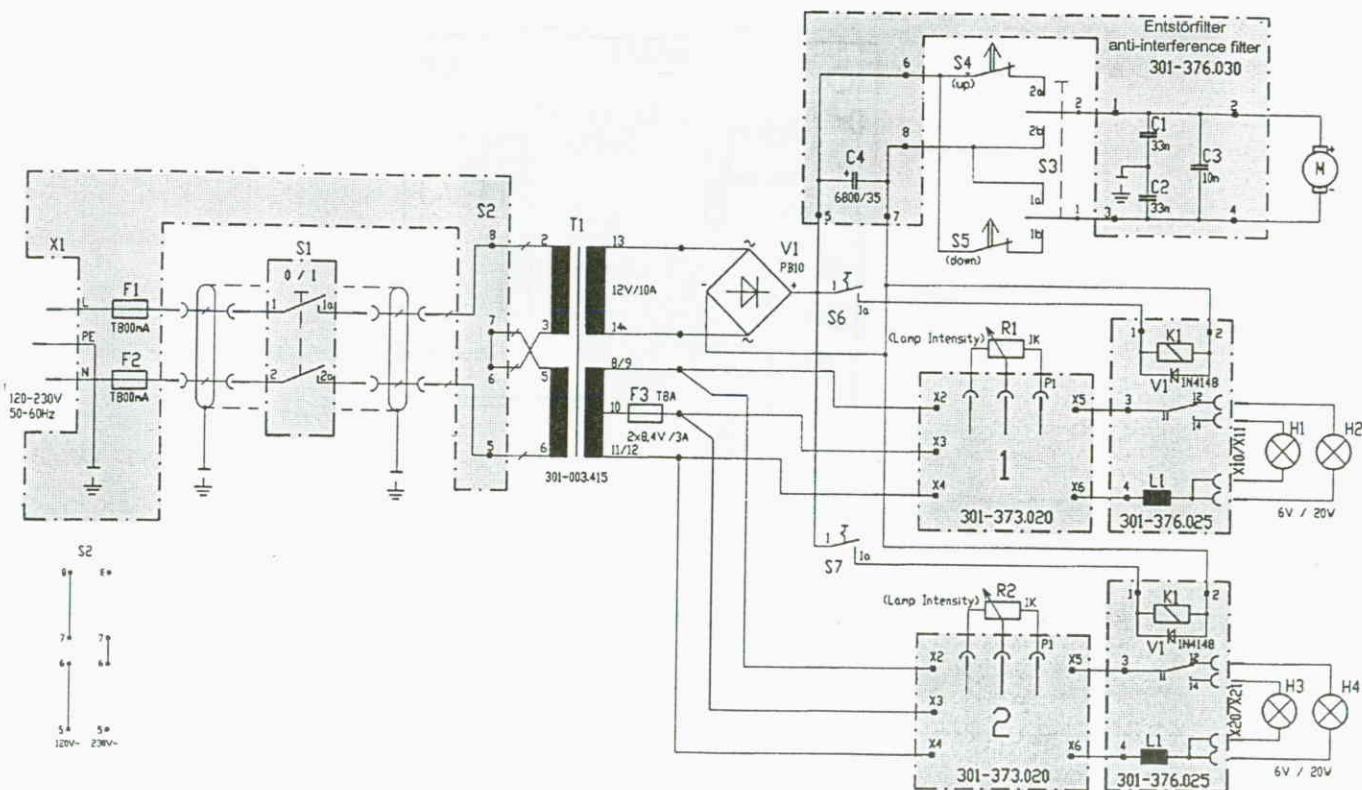
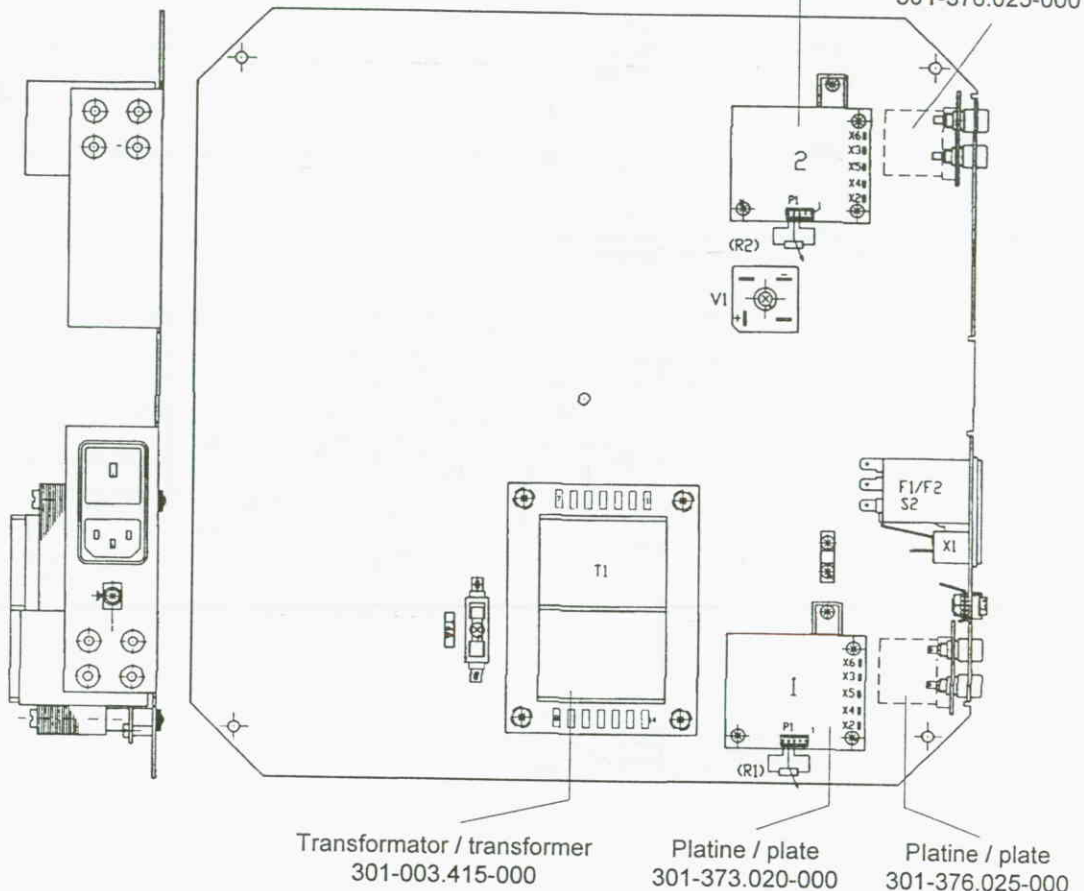
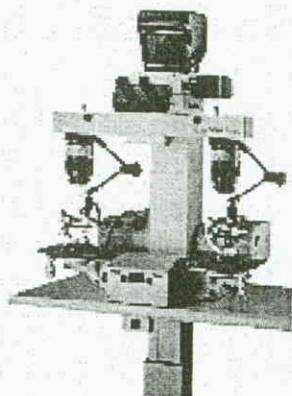
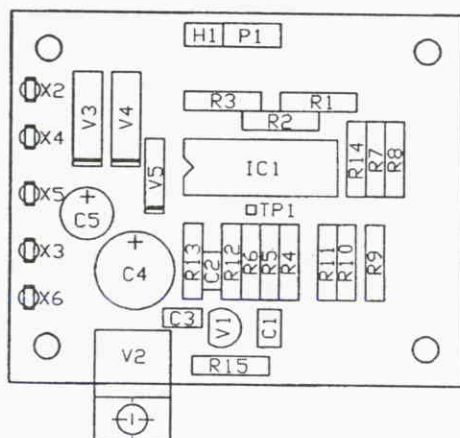
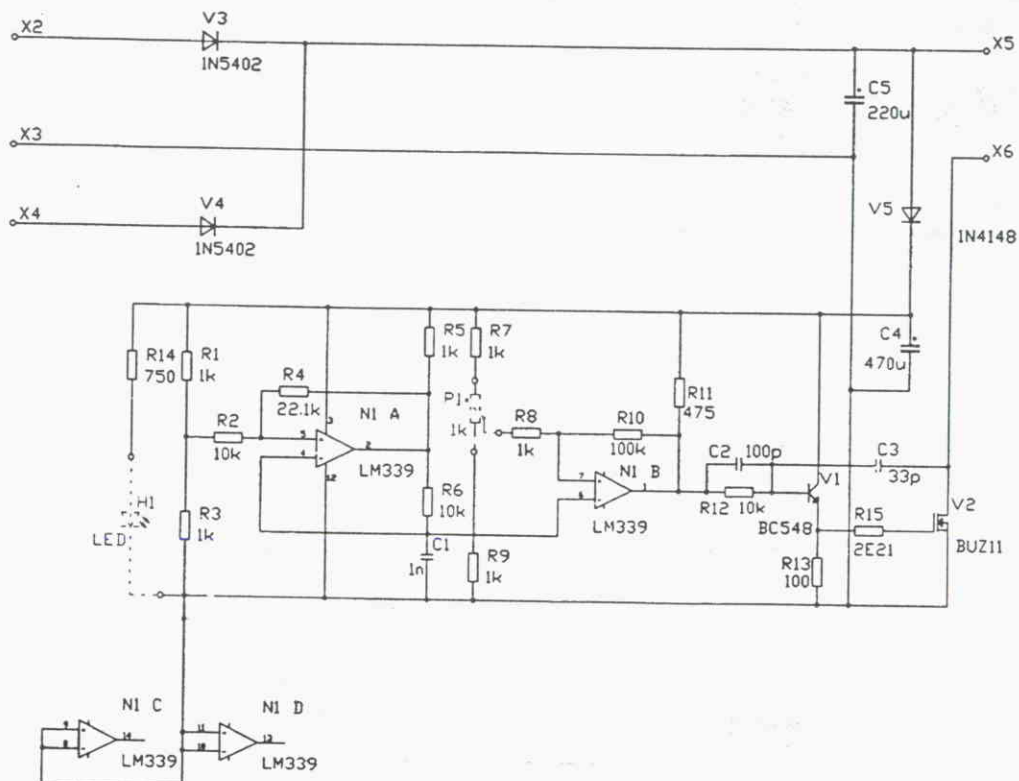
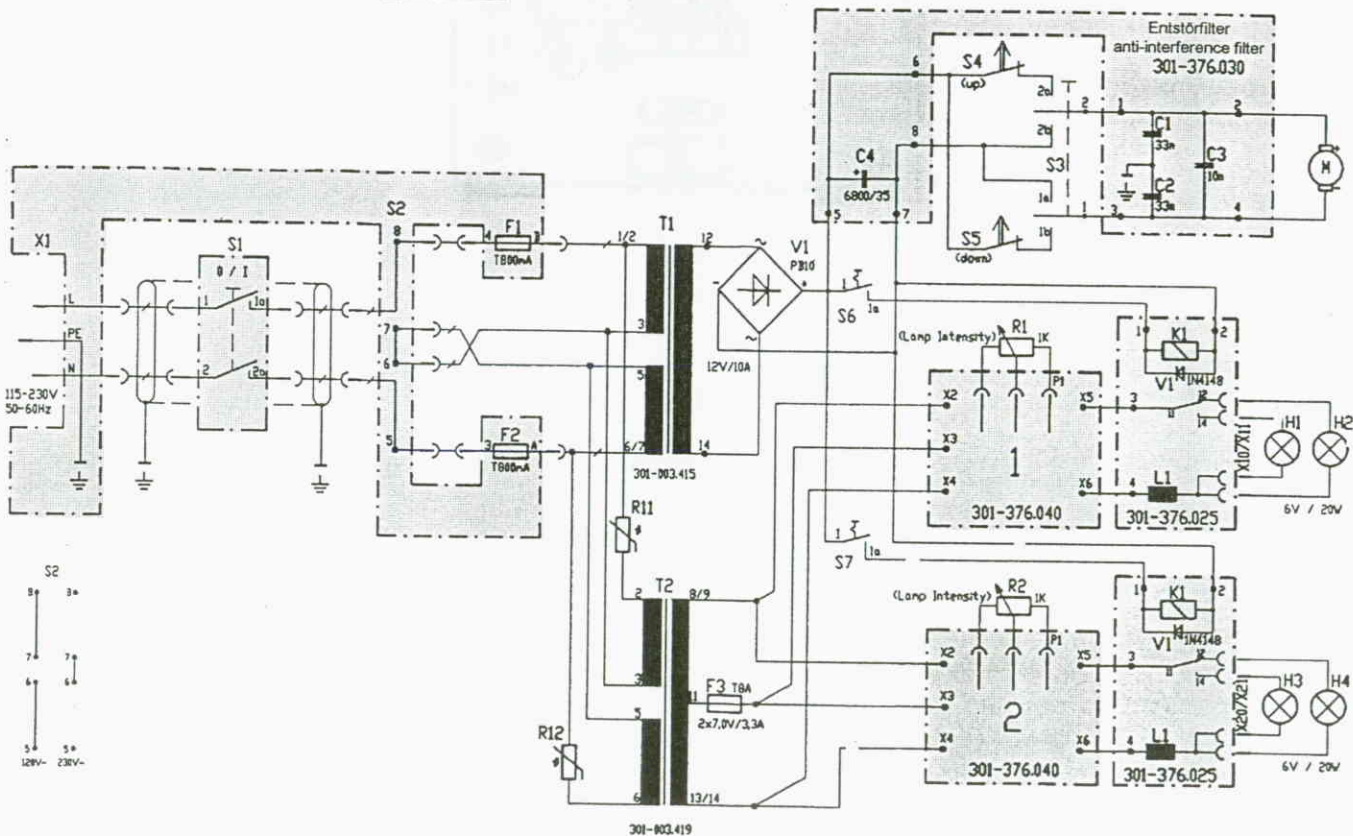
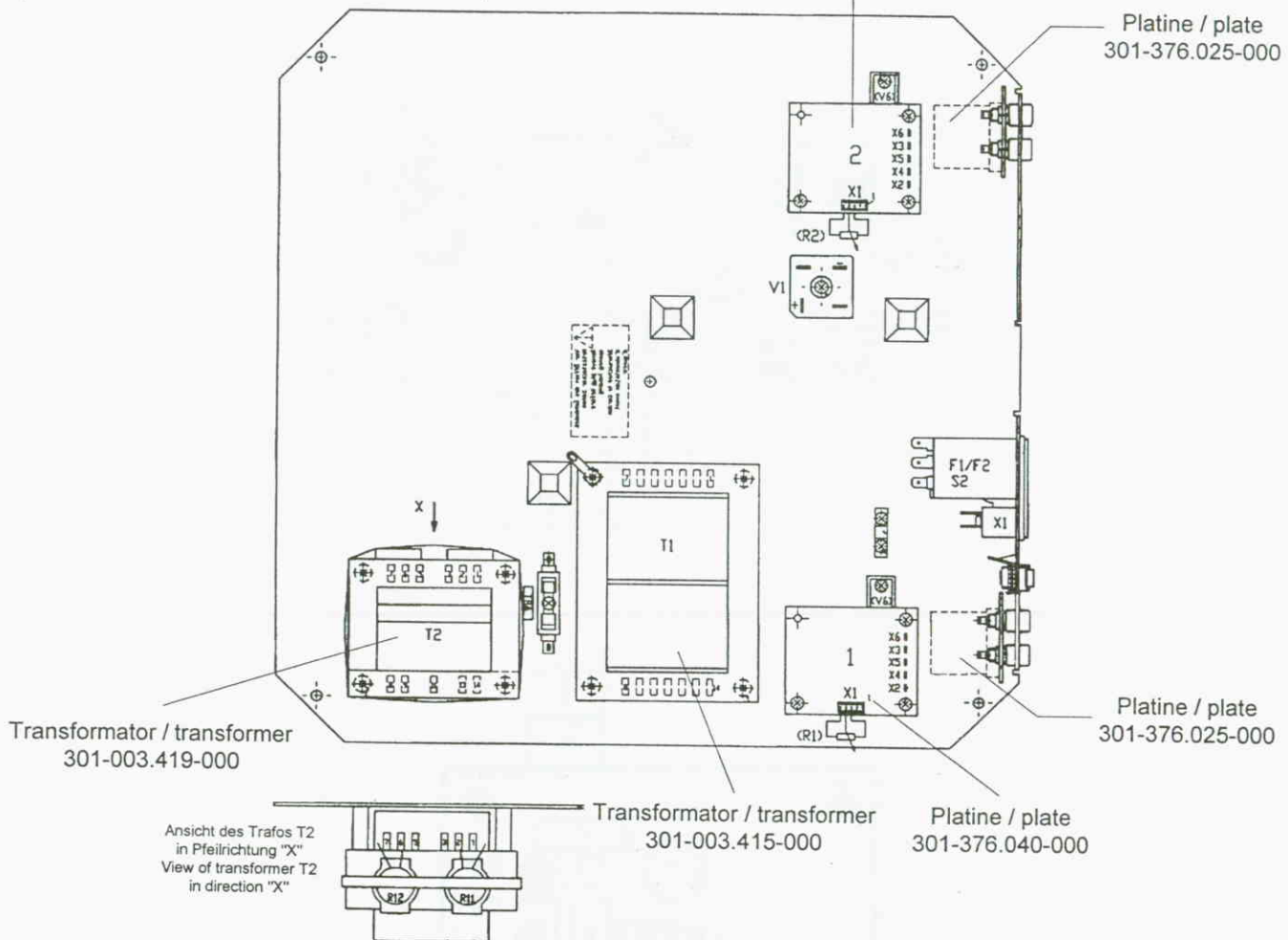
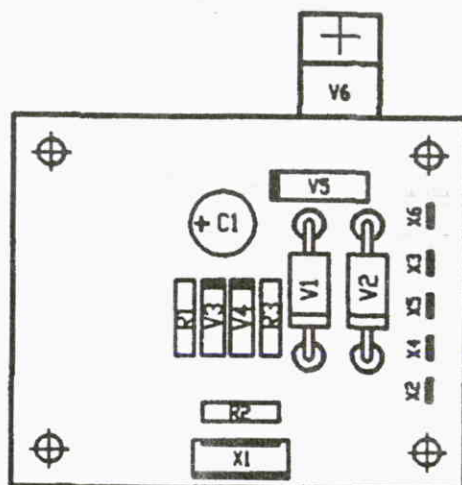
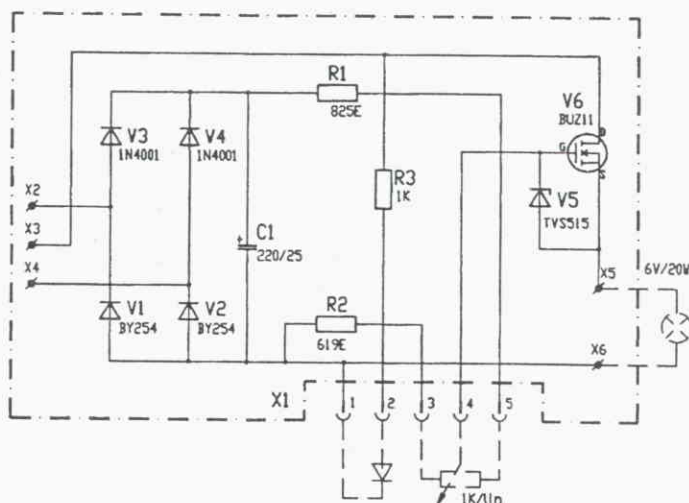
(*a) 1 Chassis Power-Supply complete 301-376.021-000

*4 Pur DC voltage (measurable with DC voltmeter)

*6 Printed circuit boards have to be exchange parallel (same manufacturer of the transistors V6 are coercive).

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Design and specifications subject to alteration without notice.

Printed in W-Germany